




Ferrule Removal Tool

- Works with PEEK™ and metal fittings
- Inexpensive and well-made
- Small enough to keep in your pocket



This tool removes high pressure PEEK™ or stainless steel ferrules that are stuck in your system. Works great for removing two-piece self-locking ferrule systems and single PEEK™ ferrules.

To use, insert the smaller drill bit end of the tool into the broken fitting. Turn clockwise to drill out a larger hole. Insert the larger end of the tool into the hole and turn counter-clockwise until a connection is felt. Gently continue turning until the fitting comes out of the port.

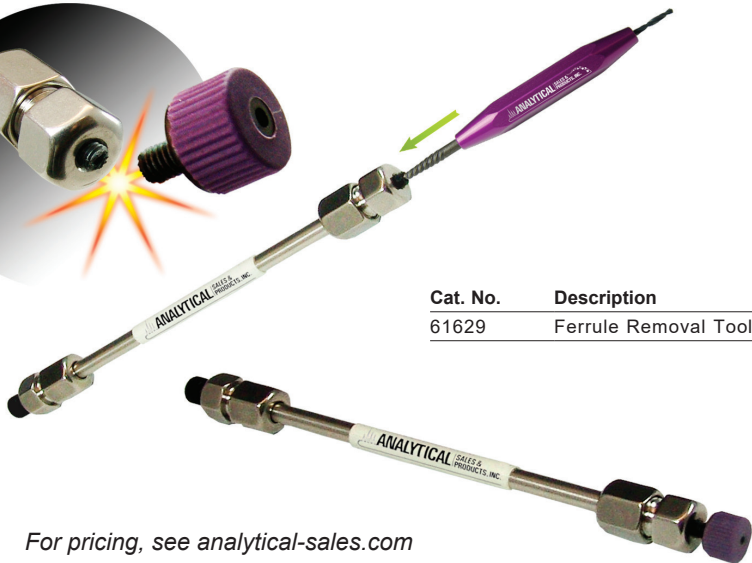

Cat. No.	Description
61629	Ferrule Removal Tool

For pricing, see analytical-sales.com