



analytical

Ferrule Removal Tool

2024

- Works with PEEK™ and metal fittings
- Inexpensive and well-made
- Small enough to keep in your pocket



This tool removes high pressure PEEK™ or stainless steel ferrules that are stuck in your system. Works great for removing two-piece self-locking ferrule systems and single PEEK™ ferrules.

To use, insert the smaller drill bit end of the tool into the broken fitting. Turn clockwise to drill out a larger hole. Insert the larger end of the tool into the hole and turn counter-clockwise until a connection is felt. Gently continue turning until the fitting comes out of the port.



| Cat. No. | Description |
|----------|----------------------|
| 61629 | Ferrule Removal Tool |

For pricing, see analytical-sales.com