

Tips for Browse Mode

How to Quickly Toggle to Your Template File.

1. Click **Power** to turn on your unit.

2. Connect the flash drive with your template files. The "USB Connected" Light will turn on.

3. Click **Mode** to toggle to "Browse" Mode.

4. Using the conversion chart below, type in the template file number with the index buttons. This will take you directly to the template file.

First Digit:	Second Digit:
A=0	1=0
B=1	2=1
C=2	3=2
D=3	4=3
E=4	5=4
F=5	6=5
G=6	7=6
H=7	8=7
I=8	9=8
J=9	10=9

Example: For template file #50, press "F" then "1".

5. Preview this template file by clicking



Or go to the next file by clicking



6. Begin running the template file by clicking to toggle to "Start Selection" Mode.



7. Step through the template file by clicking or by using the foot pedal.

