
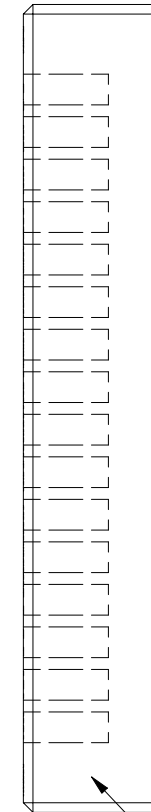
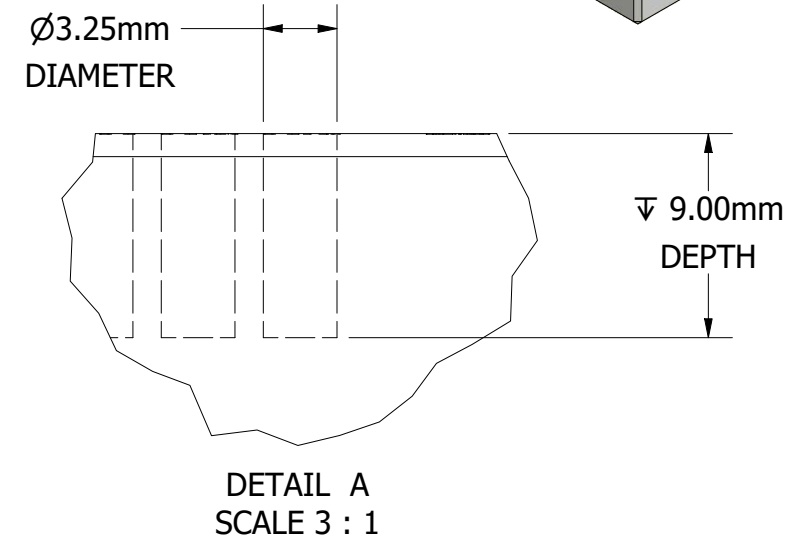
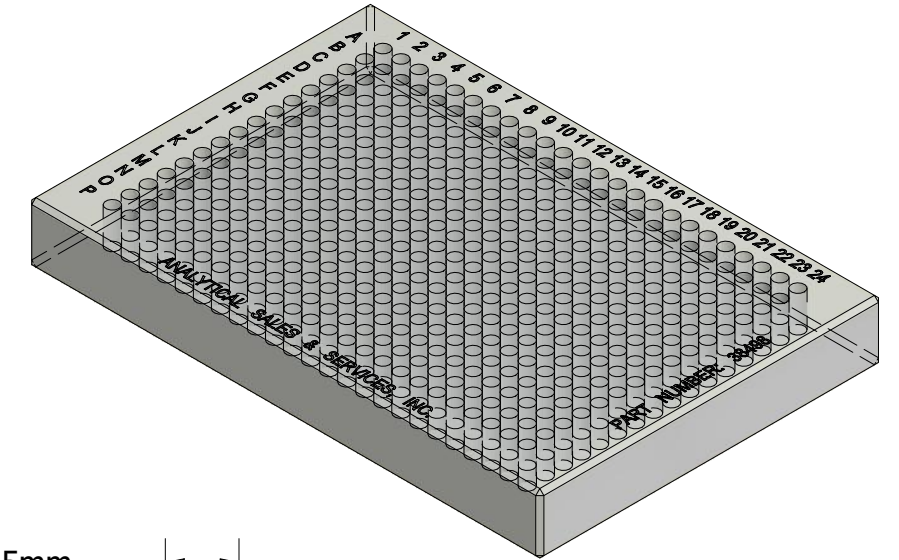


NOTES:

1. MATERIAL IS BOROSILICATE GLASS
2. PLATE POLISHED CLEAR
3. FULL CAPACITY OF WELL = 74.7µL


REVISION HISTORY			
REV	DESCRIPTION	DATE	APPROVED
1A	CUSTOMER-RELEASE PRINT	10/12/2020	ENGINEERING
1B	MADE WELL WIDER/SHALLOWER	11/5/2020	ENGINEERING
1C	ENGINEERING DIMENSIONAL ADJUSTMENT	11/30/2020	ENGINEERING

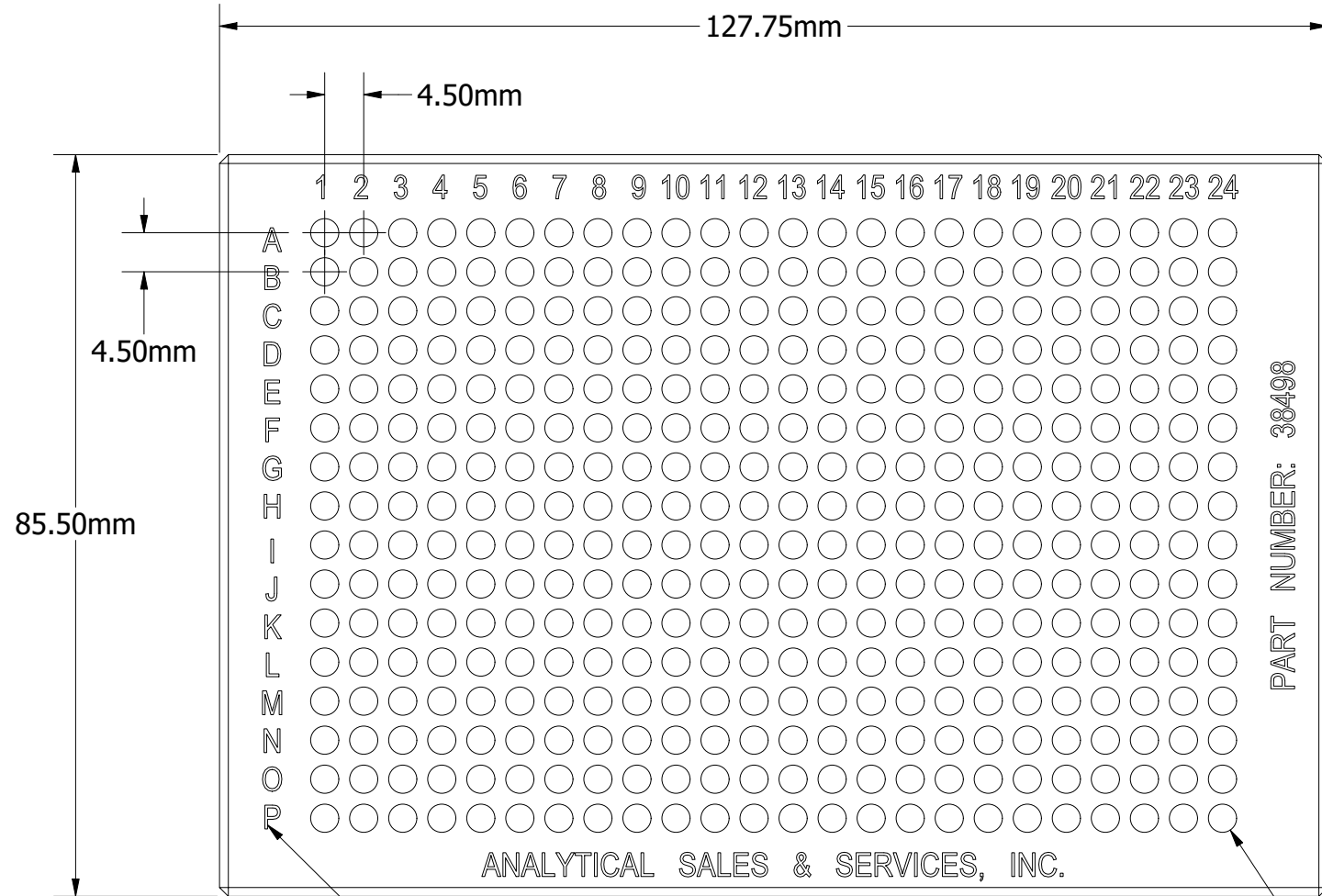

Ø3.25mm DIAMETER  
 ▽ 9.00mm DEPTH  
 (384 PLACES)

SIDE EDGES CLEAR

ALL TEXT ETCHED ONTO TOP SURFACE

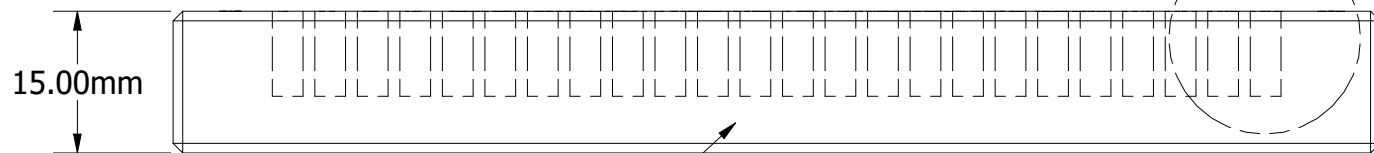

SIDE EDGES CLEAR

DRAWN Matt Dabrowski email engineering@analytical-sales.com	10/12/2020	<b>ANALYTICAL</b>	
		SALES & SERVICES, INC © 2020	
		TITLE	
		38498_384_well_glass_plate_rev1c_P	
		SIZE B	REV 1C
		SCALE 5 : 4	SHEET 1 OF 1

P

A