

Instructions for adding the EI Baidouri Shield to a DBS Punch System

1. Starting with an assembled DBS unit, loosen the punch pod from the stand rod by turning the orange handle in the punch pod clamp counterclockwise. Support the punch pod with your other hand to prevent the pod from sliding down the stand rod.
2. Slide the punch pod up to remove, and place on a clean, flat surface.
3. Once the punch pod is removed, slide the provided EI Baidouri Shield clamp assembly over the stand rod and position it so that the surface with the cutout is on top and facing the front.
4. Turn the orange handle clockwise to fasten the EI Baidouri Shield's clamp to the stand rod.
5. Position the clear EI Baidouri Shield plate so that the smaller hole matches up with the clamp's screw hole and the shield is flush with the upper surface of the clamp.
6. Insert and tighten the provided stainless steel thumbscrew until the clear EI Baidouri shield plate is securely fastened to the clamp.
7. Slide the provided silicone gasket over the glass pipette of the DBS punch pod.
8. Slide the punch pod clamp over the stand rod.
9. Position the punch pod so that the glass pipette goes through the larger hole in the EI Baidouri Shield and the silicone gasket is flush with the EI Baidouri Shield's top surface.
10. Once the punch pod is properly aligned, secure it to the stand rod by turning the orange handle on the clamp clockwise.

