



ViewOne™ LabLite / Indexer Instructions

1. Plug the LabLite/Indexer into the receptacle on the universal AC adapter (Model: 3A-061WP05, output=5V, 1.2A, input 100-240V, ~50-60Hz).
2. Plug the adapter into the power source.
3. Press the “Power” button once to turn on the LabLite and toggle to adjust intensity (high, medium, low).
4. Pressing the button while on low intensity will turn off the LabLite.

Note:

- The universal AC adapter comes with 4 different plug configurations. Before using, you must install the appropriate plug corresponding to the outlet of your country.
- Never attempt to open a ViewOne LabLite: This will compromise its function and will void any warranty.

I want another one!!!

Call, fax or email to tell us how many you want...



Catalog No.
56411

Description
ViewOne Indexer

Price
\$243



179 Route 206, Flanders, NJ 07836
P: 973-616-0700 F: 973-616-0133
www.analytical-sales.com