

**Parallel Synthesis** 

**Photoredox Catalysis** 


**Electrochemistry** 




## **Table of Contents**

## Photoredox Catalysis & Parallel Synthesis

Aluminum Reaction Blocks - Overview	2
Temperature Controlled Reactors (TCR)	4
Aluminum Reaction Blocks Micro (50µL, 4x21mm Glass Inserts): 6 • Standard (1mL, 8x30mm Vials): 8 • HPLC (2mL, 12x32mm Vials): 12 1 DRAM (4mL, 15x45mm Vials): 14 • 2 DRAM (8mL, 17x60mm Vials): 16	6
Lumidox Gen II LED arrays for Photoredox	8
Thermal Transfer Decks (TTD)2	25
Cell Culture / PCR Adapters2	26
Flow Reactor for Photoredox2	28
LumLamp for Photoredox2	29

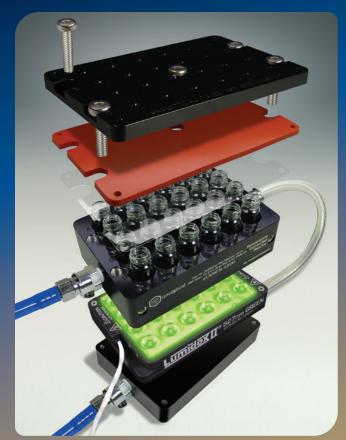



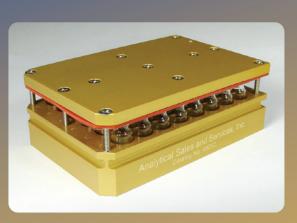





30

### **Electrochemistry**

HTe-Chem	30
Flow Flectrolysis	22

#### Accessories

EquaVAP® Evaporators: **34** • Collection Plates and Cap Mats: **36** • Thermal Adapter Plates: **37** Filter Plates and Vacuum Manifold: **38** • Vial Trays and Loaders: **39** • Powder Transfer Plates: **40** 



#### Para-dox® Aluminum Reaction Blocks

For Photoredox Catalysis and Parallel Synthesis Applications

A successful approach to high-throughput reaction screening requires the best tools available. Our comprehensive line of Photoredox Catalysis and Parallel Synthesis reaction blocks are exactly what you need. Designed specifically for high-throughput screening (HTS) applications, the SBS (SLAS/ANSI) format allows for use in any industry standard automation application.

- Useful for generating compound arrays (medicinal chemistry)
- Useful for conducting screening reaction conditions
- Can be used on tumble stirrers / hot plate stirrers / robotic platforms
- Can easily be used in a glovebox
- Temperature range -78°C 150°C
- Validated to have less than 5% solvent loss with prolonged heating above boiling point
- Silicone Rubber Mats provide compression sealing
- Teflon® PFA Films keep the glass reaction vials from sticking to the silicone rubber mats during heating

#### For 4x21mm (50µL) Flat Bottom Glass Inserts

Gen II







Micro 96-Position Photoredox Catalysis, Gen II Photoredox Catalysis, Gen II



Micro 24-Position **Photoredox Catalysis** 



Pages 6, 7

**Pages 8-11** 

**Micro 96-Position Photoredox Catalysis** 

#### For 8x30mm (1mL) Clear Glass Shell or Crimp Top Vials



Standard 96-Position Photoredox Catalysis, Gen II



Standard 96-Position Parallel Synthesis, Gen II



Standard 96-Position

24249



Standard 96-Position Photoredox LIGHTWEIGHT Photoredox/Parallel Synthesis



84001-Case<sup>3</sup>

10421-Case\*

Gen II

**Standard 24-Position Photoredox** 



Catalysis/Parallel Synthesis, Gen II



Standard 24-Position **Photoredox Catalysis/Parallel Synthesis** 



48-Position TCR

#### For 12x32mm (2mL) Crimp/Snap Vials




Photoredox Catalysis, Gen II



**HPLC 48-Position** Parallel Synthesis, Gen II



**HPLC 48-Position Photoredox Catalysis** 

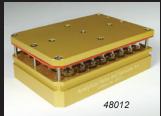

Pages 12, 13

Pages 14, 15

**HPLC 48-Position Parallel Synthesis** 

#### For 15x45mm (4mL) Clear Glass Screw Vials



1 Dram 24-Position Photoredox Catalysis, Gen II (18mm well spacing)

1 Dram 24-Position

**Photoredox Catalysis** 


(20mm well spacing)




1 Dram 24-Position Parallel Synthesis, Gen II (18mm well spacing)



1 Dram 24-Position Photoredox Catalysis, Gen II w/ Temperature Transfer Cover

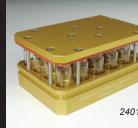

1 Dram 24-Position **Parallel Synthesis** (20mm well spacing)

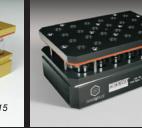

1 Dram 24-Position **Photoredox Catalysis** (18mm well spacing)



24-Position TCR

#### For 17x60mm (8mL) Clear Glass Screw Vials

31531-Case



2 Dram 24-Position Photoredox Catalysis, Gen II (18mm well spacing)



2 Dram 24-Position Parallel Synthesis, Gen II (18mm well spacing)



2 Dram 24-Position **Photoredox Catalysis** 



Pages 16, 17

2 Dram 24-Position **Parallel Synthesis** 

www.analytical-sales.com \* See page 39 for Vial Loader Trays


31760-Case

#### **Photoredox Catalysis** & Parallel Synthesis




#### **Temperature Controlled Reactors (TCR) for** 8x30mm (1mL) and 15x45mm (1 Dram) Vials

- · Provides extremely uniform thermal control for high throughput experimentation (HTE) with a temperature difference of +/- 1°C
- Capable of enhanced reproducibility in heating and cooling applications
- High quality leak-proof CPC fittings for quick and easy tubing connection and removal
- Compatible with a wide range of heat transfer fluids such as water (down to 5°C), ethylene glycol, polypropylene glycol and siliconebased fluids (ie SYLTHERM™)
- Designed to standard SLAS dimensions (127.75mm x 85.5mm)
- Compatible with auto-samplers and high throughput chemistry devices
- 4mm holes in lid allow for use with most common autosampler needles
- Threaded holes (6-32) in lid for accessory attachment

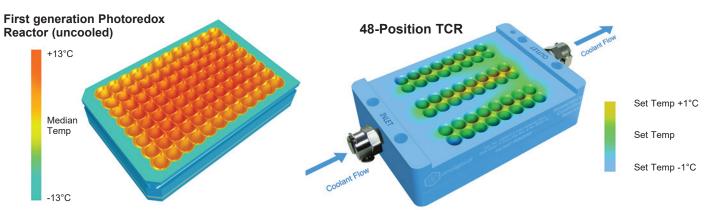

Simulated heat maps of a standard 96-position Photoredox Reactor Block vs. a 48-position Temperature Controlled Reactor

(TCR) when used with a Lumidox®II 48 Position LED Array at full power (stage 5).

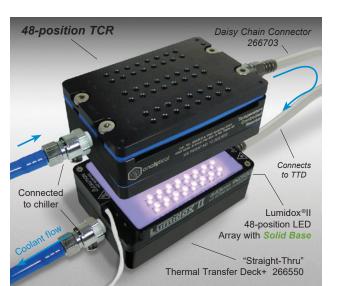

Typical configuration: 48-position TCR used in conjunction with a Lumidox®II 48-position Solid Base LED Array, a "Straight-Thru" Thermal Transfer Deck+ (for additional cooling of array) and connected to an external liquid chiller

#### **48-Position Temperature Controlled Reactor**

• For 8x30mm (1mL) Vials • 48 Wells (6 Rows of 8) • 9mm Well Spacing


Cat. No.	Description	Qty
489900	48-position TCR. Includes: PFA Film, Blue Fluorosilicone Sealing Mat,	Each
	Rubber Gasket, Silicone O-Rings, CPC Fittings, PUR Tubing and Screws.	
489906	1/8" Thick, Blue Fluorosilicone Rubber Sealing Mats for TCR	5
489907	0.005" Thick PFA Sealing Films for TCR	25
489908	9mm OD, 6mm ID, 1.5mm Wide Silicone O-rings for TCR	50
SHS0034	SS Low-Profile Socket Head Screw w/ Hex Drive, 5/16"-18 Threads, 3/4" Long	5
84001-Case	1mL Clear Glass Shell Vials, 8 x 30mm	1000
488401	Vial Tray for TCR, Pre-loaded with 48 8x30 Shell Vials (84001-CASE)	Ea

#### Solid Base Lumidox®II 48-position LED Array for TCR

- Lens Mat surface
- Requires Thermal Transfer Deck (TTD) connected to chiller

Cat. No.	Wavelength	Cat. No.	Wavelength
LUM248LS365	UV365	LUM248LS405	UV405
LUM248LS375	UV375	LUM248LS445	445-INDIGO
LUM248LS385	UV385	LUM248LS470	470-BLUE
LUM248LS395	UV395	LUM248LS505	505-CYAN

1	Cat. No.	Wavelength
5	LUM248LS527	527-GREEN
)	LUM248LS590	590-AMBER
	LUM248LS630	630-RED
1	LUM248LSWHT	WHITE

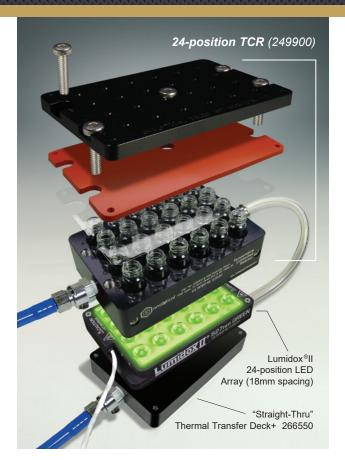



#### Screen & Scale-up with the Same Precision!

#### 24-Position Temperature Controlled Reactor

- For 15x45mm (1 Dram) Vials
- 24 Wells (4 Rows of 6)
- 18mm Well Spacing

at. No.	Description	Qty
49900	24-position TCR. Includes: PFA Film, Silicone Rubber Sealing Mat,	Each
	Rubber Gasket, CPC Fittings, PUR Tubing, Screws. Vials not included.	
01247	Top Rubber Mats for Gen II 24-Position Reactors (with 18mm Spacing)	25
01248	Top PFA Films for Gen II 24-Position Reactors (with 18mm Spacing)	25
SHS0114	Socket Head Screws for Gen II 24-Position Reactors, 5/16"-18 Threads, 1-1/4" Long	5
1531-Case	13mm, 15 x 45mm, 4mL (1 Dram) Clear Glass Screw Vials	1000
1554-Case	13mm Black PP Solid Screw Caps with PTFE/F217 Liners	1000

#### Lumidox®II 24-position LED Array with 18mm Spacing, Solid Base

· Lens Mat surface

LU

• Requires Thermal Transfer Deck (TTD) connected to chiller

at. No.	Wavelength (nm)	Cat. No.	Wavelength (nm)
UM22418LS255	UVC-255	LUM22418LS445	445-INDIGO
UM22418LS365	UV365	LUM22418LS470	470-BLUE
UM22418LS375	UV375	LUM22418LS505	505-CYAN
UM22418LS385	UV385	LUM22418LS527	527-GREEN
UM22418LS395	UV395	LUM22418LS590	590-AMBER
UM22418LS405	UV405	LUM22418LS630	630-RED
UM22418LS420	420-VIOLET	LUM22418LSWHT	WHITE

#### 24-Position TCR with Temperature Transfer Cover

Perfect for those who prefer to cap their vials instead of sealing them with the standard aluminum lid, rubber mat and PFA film that come with the block assembly. When used in conjunction with a Thermal Transfer Deck+ (TTD+) and external liquid chiller, the cover helps control excessive heat generated by a Lumidox® II LED array and also greatly reduces light bleed.

The **Temperature Transfer Cover** is also available separately (249913) and can be used with other Para-dox® Gen II 24-position reactor blocks (18mm spacing) for photoredox and parallel synthesis.

Cat. No.	Description	Qty
249800	24-Position TCR with Temperature Transfer Cover	Each
249913	Temperature Transfer Cover for Gen II, 24-Position, 18mm Spacing Reactors	Each

#### **Additional Required Parts**


Cat. No.	Description	Qty
266550	Para-dox <sup>®</sup> "Straight-Thru" Thermal Transfer Deck (TTD)	Each
266703	Daisy Chain Connector for TCR and "Straight-Thru" TTD	Each
TCubeEdge	Recirculating Chiller. Operating Range: 0°C - 65°C	Each



#### Tech Tip

- Temperature Controlled Reactor (TCR)\* cools vials
- Thermal Transfer Deck (TTD)\* cools LED Array (solid base)
- LED Array with solid base needs TTD and chiller for cooling
- \* Connects to External Liquid Chiller, required



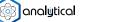

## Photoredox Catalysis & Parallel Synthesis



## Micro 24-Position Photoredox Catalysis Aluminum Reaction Blocks, Gen II

For 4x21mm, 50µL Glass Inserts

Catalog No.	Description	Qty
107240	Gen II, 24 Position, 4-BOLT, Photoredox Reaction Block for 4x21mm (50μL) Vials, with 9mm Well Spacing	Each
107245	Bottom Rubber Mats for 107240, Gen II 24 Position Reactor (4x21mm vials, 9mm spacing)	25pk
102247	Top Rubber Mats for Gen II 24 Position Reactor (9mm spacing)	25pk
102248	Top PFA Films for Gen II 24 Position Reactor (9mm spacing)	25pk
SHS0058	10-24, 5/8" Long, SS Socket Head Screw	5







#### Micro 96-Position Photoredox Catalysis Aluminum Reaction Blocks, Gen II For 4x21mm, 50µL Glass Inserts

Catalog No.	Description	Qty
103960	Gen II, 96 Position, 4-BOLT, Photoredox Reaction Block for 4x21mm (50μL) Vials, with 9mm Well Spacing	Each
103965	Bottom Rubber Mat for 103960, Gen II 96 Position Reactor (4x21mm vials, 9mm spacing)	Each
101967	Top Rubber Mats for Gen II 96-Position and 48-Position Reactors	25
101968	Top PFA Films for Gen II 96-Position and 48-Position Reactors	25
SHS0034	Socket Head Screws for Gen II 96-Position and 48-Position Reactors	5







## TECH NOTE: Replace PFA Films with every use for best performance

## Gen II Gen II NEW! Para-dox® Gen II Aluminum Reaction Blocks

#### For Photoredox Catalysis and Parallel Synthesis

#### Benefits of the Gen II line include:

- Less bolts = less time to assemble
- Less light bleed when used with Lumidox<sup>®</sup> II LED arrays (48 and 96 position blocks)
- Less Maintenance

Gen II

- Vials fit tighter (less shifting)
- Longer lasting screws (larger diameter, larger threads)

#### Additional features:

- · Port for thermocouple, sensors, etc.
- Corner holes on both the lid and the bottom for accessory attachment
- Larger holes in lid to accommodate a variety of needle sizes
- Uses only one 1/8" rubber mat compared to two 1/16" mats on previous versions (makes for easier assembly)










## Micro 24-Position Photoredox Catalysis Aluminum Reaction Blocks

For 4x21mm, 50µL Glass Inserts

Cat. No.	Description	Qty
24250	Photoredox 24-Position Micro Block Assembly. Includes Vial Rack, Bottom and Top Covers, PFA Sheets, Rubber Mats and Screws	Each
24257	Bottom PFA Films	25
24258	Bottom Rubber Mats	25
24261	Top PFA Films for 24-Position Photoredox Micro	25
24262	Top Rubber Mats for 24-Position Photoredox Micro	25
1/2"SCREW96	1/2" Bottom Screws	100
1"SCREW	1" Bottom Screws	100

 4x21mm flat-bottom glass micro vials, 50µL volume (25µL maximum volume if using tumble stir bars)



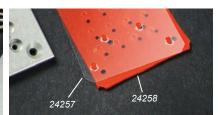

#### Micro 96-Position Photoredox Catalysis Aluminum Reaction Blocks

For 4x21mm, 50µL Glass Inserts

Description	Qt
Photoredox 96-Position Micro Block Assembly. Includes Vial Rack,	Eac
Bottom and Top Covers, PFA Sheets, Rubber Mats and Screws	
Bottom PFA Films	2
Bottom Rubber Mats	2
Top PFA Films for 96-Position Photoredox Micro	2
Top Rubber Mats for 96-Position Photoredox Micro	2
1/2" Bottom Screws	10
1" Top Screws	10
	Photoredox 96-Position Micro Block Assembly. Includes Vial Rack, Bottom and Top Covers, PFA Sheets, Rubber Mats and Screws Bottom PFA Films Bottom Rubber Mats Top PFA Films for 96-Position Photoredox Micro Top Rubber Mats for 96-Position Photoredox Micro 1/2" Bottom Screws










www.analytical-sales.com

#### Accessories


Cat. No.	Description	Qty
10421-Case	Micro Vials - 50µL Flat Bottom Glass Inserts, 4x21mm	1000
13257	Magnetic Tumble Stir Bars, 1.32mm x 1.57mm (Fleas)	1000
SD1000	Milwaukee M4 1/4in. Hex Screwdriver Kit. Includes 2 Batteries and Charger (Note: batteries not available for international shipping)	Each

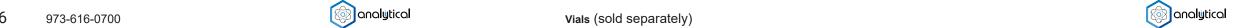

#### **Photoredox Catalysis** & Parallel Synthesis





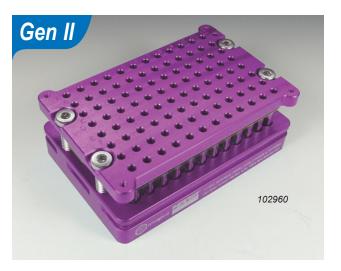


#### 96-Position Reactor Assembly Weight\* Gen II, Lightweight (102960): 477.8g - Base only, no lid or screws: 267.8g 769.0g Gen II, 4-Bolt (101960): Gen I, Original (96973): 536.0g

\* Weight without vials. Adding vials will add ~ 131g.

973-616-0700



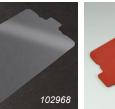



analytical





#### **Standard 96-Position Photoredox Catalysis Aluminum Reaction Blocks, Gen II**


for 1mL, 8x30mm Glass Inserts

Catalog No.	Description	Qty
101960	Gen II, 96 Position, 4-BOLT, Photoredox Reaction Block for 8x30mm (1mL) Vials, with 9mm Well Spacing	Each
101965	Bottom Rubber Mats for Gen II 96-Position Reactor (101960)	25
101964	Bottom PFA Film for Gen II 96-Position Reactor (101960)	Each

#### **Standard 96-Position Parallel Synthesis** Aluminum Reaction Blocks, Gen II

for 1mL, 8x30mm Glass Inserts

Catalog No.	Description	Qty
104960	Gen II, 96 Position, 4-BOLT, Parallel Synthesis Reaction Block for 8x30mm (1mL) Vials, with 9mm Well Spacing	Each
104965	Bottom Rubber Mats for Gen II 96 Position Reactor (104960)	25PK
104964	Bottom PFA Film for Gen II 96-Position Reactor (104960)	Each

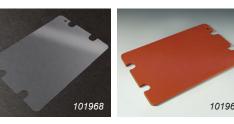



#### Replacement Parts (for photoredox AND parallel synthesis Gen II 96-position blocks)

Catalog No.	Description	Qty
101967	Top Rubber Mats for Gen II 96-Position and 48-Position Reactors	25
101968	Top PFA Films for Gen II 96-Position and 48-Position Reactors	25
SHS0034	Socket Head Screws for Gen II 96-Position and 48-Position Reactors	5

#### Standard 96-Position Photoredox LIGHTWEIGHT Reaction Block, Gen II for 1mL, 8x30mm Glass Inserts

- 37% lighter than Gen II, 96-position standard 4-bolt reactor
- Great for use on a low-capacity balance with a 500g max limit
- Reduced weight allows for better centrifuge performance
- 9mm Well Spacing

Catalog No.	Description	QTY
102960	Gen II, 4-Bolt, 96-Position LIGHTWEIGHT Photoredox Reaction Block for 1mL, 8x30mm Glass Inserts. Includes Vial Rack, Top Cover, PFA Sheet, Rubber Mats and Screws	Each
102968	Top PFA Films for Gen II LIGHTWEIGHT 96-Position Reactor (102960)	25
102967	Top Rubber Mats for Gen II LIGHTWEIGHT 96-Position Reactor (102960)	25
102965	Bottom Rubber Mat for Gen II LIGHTWEIGHT 96-Position Reactor (102960) - (two required per block)	Each
SHS0114	Socket Head Screws for Gen II 24-Position Reactors and 96-Position Lightweight Reactors	5

## Gen



#### **Standard 96-Position Photoredox Catalysis Aluminum Reaction Blocks**


for 1mL, 8x30mm Glass Inserts

at. No.	Description	Qty
6973	Photoredox 96-Position Block Assembly Includes: Vial Rack, Covers, Mats, Films and Screws	Each
6979	Replacement Bottom Films for 96-Position Photoredox Block	25
6980	Replacement Bottom Mats for 96-Position Photoredox Block	25

#### **Standard 96-Position Parallel Synthesis Aluminum Reaction Blocks**

for 1mL, 8x30mm Glass Inserts

Cat. No.	Description	Qty
96960	Parallel Synthesis 96-Position Block Assembly. Includes Vial Rack, Top and Bottom Covers, PFA Films, Rubber Mats and Screws	Each

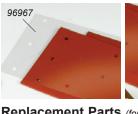





#### Replacement Parts (for photoredox AND parallel synthesis 96-position blocks)

Cat. No.	Description	Qty
6967	Replacement Films for 96-Position Optimization Block	25
6965	Replacement Mats for 96-Position Optimization Block	25
/2"SCREW96	1/2" Bottom Screws	100
"SCREW	1" Top Screws	100

#### Accessories

Clear, Round Well Cap Mat with Screw Holes (allows for assembly with catalysis reaction blocks)

Cat. No.	Description	Qty
964007H	Clear Ultra Thin Round Well Cap Mat,	5
	Silicone/PTFE, with Screw Holes	

#### Plugs for 8x30mm Vials Inserts

Cat. No.	Description	Qty
96077	Pierceable TPE Plugs on 96 Well Spine,	5
	Individual, for Tubes/Vials, Purple	
96089	Pierceable TPE Plugs on 96 Well Spine,	5
	Strips of 8, for Tubes/Vials, Natural	




#### **Top Cover with Wider Holes for 96-Position Reaction Blocks**

- Enlarged (4.3mm) Holes
- 316 Stainless Steel
- Made for Automated Sampling Instruments

#### Cat. No. Description

96985 Reaction Block Cover w/ 4.3mm Holes, 6mm Thick 96986 Reaction Block Cover w/ 4.3mm Holes, 3mm Thick

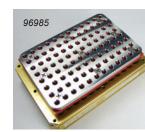

Gen

#### Barcoded 8x30mm Vial Inserts in Tray for 96-Well Blocks

Cat. No. Description 9984001 Barcoded 8x30mm Vial Inserts in Tray for 96-Well Blocks

#### Stir Bars/Power Screwdriver


Cat. No.	Description	Qty
13258	SS Cylinder Stir Bars, 1.98mm x 4.80mm (for 8x30mm Vials)	1000
SD1000	Milwaukee M4 1/4in. Battery Powered Hex Screwdriver Kit	Ea



#### Spacers for 5x31mm Inserts

Allows 250µL, 5x31mm inserts (part no. 20303) to be used in 24 and 96 position blocks for 8x30mm inserts.

Cat. No.	Description	Qty
96969	Anodized Aluminum Spacers	96
	for 5x31mm Inserts	
20303-Case	250μL, 5 x 31mm Inserts	1000

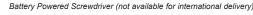

#### **Photoredox Catalysis** & Parallel Synthesis







#### **Standard 24-Position Photoredox Catalysis Aluminum Reaction Blocks, Gen II**

for 1mL, 8x30mm Glass Inserts

Catalog No.	Description	Qty
102240	Gen II, 24 Position, 4-BOLT, <b>Photoredox</b> Reaction Block for 8x30mm (1mL) Vials, with 9mm Well Spacing	Each
102245	Bottom Rubber Mats for 102240, Gen II 24 Position Reactor (8x30mm vials, 9mm spacing)	25

#### **Standard 24-Position Parallel Synthesis** Aluminum Reaction Blocks, Gen II

for 1mL, 8x30mm Glass Inserts

Catalog No.	Description	Qty
105240	Gen II, 24 Position, 4-BOLT, <b>Parallel Synthesis</b> Reaction Block for 8x30mm (1mL) Vials, with 9mm Well Spacing	Each
105245	Bottom Rubber Mats for 105240, Gen II 24 Position Reactor (8x30mm vials, 9mm spacing, Closed)	25







Replacement Parts (for photoredox AND parallel synthesis Gen II 24-position blocks)

Catalog No.	Description	Qty
102247	Top Rubber Mats for Gen II 24 Position Reactors (with 9mm Spacing)	25
102248	Top PFA Films for Gen II 24 Position Reactors (with 9mm spacing)	25
SHS0058	10-24, 5/8" Long, SS Socket Head Screw	5

#### Gen II Rack Adapter for 24 Well, 9mm Spacing Gen II Reactors

(for photoredox AND parallel synthesis blocks)

Catalog No.	Description	Qty
102249	Gen II Rack Adapter for 24 Well, 9mm Spacing Gen II Reactors	Each

#### Parallel Synthesis vs. Photoredox Catalysis **Reaction Blocks**

Reactors for both applications are nearly identical. The difference lies in the base (vial rack) of the block. Reactors for parallel synthesis have wells with closed bottoms, whereas photoredox blocks have wells with open holes in the bottom allowing for light transmission via a Lumidox®II Array, as well as to read 2D barcodes.











24240

Gen I

Gen


TECH NOTE:

Replace PFA Films with every use for best performance

#### **Standard 24-Position Photoredox Catalysis Aluminum Reaction Blocks**

for 1mL, 8x30mm Glass Inserts

Cat. No.	Description	Qty
24253	Photoredox 24-Position Block Assembly. Includes Vial Rack, Top Cover, Bottom Cover, PFA Films, Rubber Mats, Screws	Each
24259	Replacement Bottom Films for 24 Position, Photoredox Block	25
24260	Replacement Bottom Mats for 24 Position, Photoredox Block	25

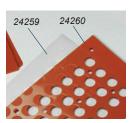


assembly (24253) only

#### **Standard 24-Position Parallel Synthesis Aluminum Reaction Blocks**


for 1mL, 8x30mm Glass Inserts

Cat. No.	Description	Qty
24240	Parallel Synthesis 24-Position Block Assembly. Includes Vial Rack, Top Cover, Bottom Cover, PFA Films, Rubber Mats, Screws and Rack Adapter	Each
24249	Parallel Synthesis/Optimization 24-Position Block Assembly, no Rack Adapter	Each

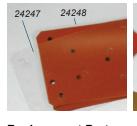






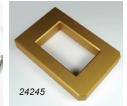

#### Replacement Partsv (for photoredox AND parallel synthesis 24-position blocks)

Cat. No.	Description	Qty
24247	Replacement Films for 24 Position, Optimization Block	25
24248	Replacement Mats for 24 Position, Optimization Block	25
3/8"SCREW24	3/8" Bottom Screws	100
1"SCREW	1" Top Screws	100
24245	Rack Adapter, Optional	Each



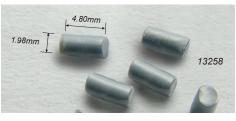


#### **Accessories**

24249

Cat. No.	Description	Qty
84001-Case	1mL Clear Glass Shell Vials, 8 x 30mm	1000
13258	Stainless Steel Cylinder Stir Bars, 1.98mm x 4.80mm (for 8x30mm Vials)	1000
SD1000	Milwaukee M4 1/4in. Hex Screwdriver Kit. Includes 2 Batteries and Charger	Each
	(Note: batteries not available for international shipping)	



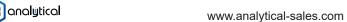

## **HPLC 48-Position Photoredox Catalysis Aluminum Reaction Blocks, Gen II**

HPLC (2mL, 12x32mm VIALS), GEN II

for 2mL, 12x32mm Vials

Catalog No.	Description	Qty
101480	Gen II, 48 Position, 4-BOLT, Photoredox Reaction Block for 12mm OD	Each
	(2mL) Vials, with 13mm Well Spacing	

## **HPLC 48-Position Parallel Synthesis Aluminum Reaction Blocks, Gen II**

for 2mL, 12x32mm Vials

Catalog No.	Description	Qty
104480	Gen II, 48 Position, 4-BOLT, Parallel Synthesis Reaction Block for 12mm OD	Each
	(2mL) Vials, with 13mm Well Spacing	









Replacement Parts (for photoredox AND parallel synthesis Gen II 48-position blocks)

Catalog No.	Description	Qty
101968	Top PFA Films for Gen II 96-Position and 48-Position Reactors	25
101967	Top Rubber Mats for Gen II 96-Position and 48-Position Reactors	25
SHS0034	Socket Head Screws for Gen II 96-Position and 48-Position Reactors	5

#### Gen II Gen II

#### **NEW!** Para-dox® Gen II Aluminum Reaction Blocks

#### For Photoredox Catalysis and Parallel Synthesis

#### Benefits of the Gen II line include:

- Less bolts = less time to assemble
- Less light bleed when used with Lumidox<sup>®</sup> II LED arrays (48 and 96 position blocks)
- Less Maintenance
- · Vials fit tighter (less shifting)
- Longer lasting screws (larger diameter, larger threads)

#### Additional features:

- Port for thermocouple, sensors, etc.
- Corner holes on both the lid and the bottom for accessory attachment
- · Larger holes in lid to accommodate a variety of needle sizes
- Uses only one 1/8" rubber mat compared to two 1/16" mats on previous versions (makes for easier assembly)

### Parallel Synthesis vs. Photoredox Catalysis Reaction Blocks


Reactors for both applications are nearly identical. The difference lies in the base (vial rack) of the block. Reactors for parallel synthesis have wells with closed bottoms, whereas photoredox blocks have wells with open holes in the bottom allowing for light transmission via a Lumidox®II Array, as well as to read 2D barcodes.

#### Parallel Synthesis Block, Gen II



Photoredox Catalysis Block, Gen II



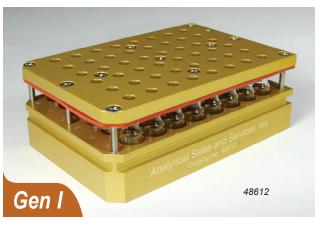

## Gen I 48012

## **HPLC 48-Position Photoredox Catalysis Aluminum Reaction Blocks**


for 2mL, 12x32mm Vials


Catalog No.	Description	Qty
48612	48-Position <b>Open Top/Open Bottom</b> Aluminum Reaction Block for 12mm (2mL) OD Vials. Includes: Base Plate, Cover, Screws, PFA Film and 2 Rubber Mats	Each

## HPLC 48-Position Parallel Synthesis Aluminum Reaction Blocks

for 2mL. 12x32mm Vials

Catalog No.	Description	Qty
48012	48-Position Aluminum Reaction Block for 12mm (2mL) OD Vials. Includes: Base Plate, Cover, Screws, PFA Film and 2 Rubber Mats	Each

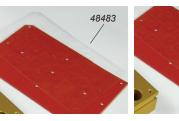

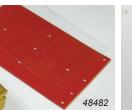




Replacement Parts (for photoredox AND parallel synthesis 48-position blocks)

Catalog No.	Description	Qty
48482	Rubber Mats for 48 Position (12mm OD) Reaction Blocks	25
48483	PFA films for 48 Position (12mm OD) Reaction Blocks	25
VSCREW48	1 1/4" Screws for 48-Position Aluminum Reaction Plate	100







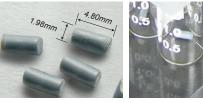

#### SS Cylinder Stir Bars for 2mL HPLC Vials

- Parylene Encapsulated
- For Photoredox and Parallel Synthesis Reaction Block Systems

Cat. No.	Description	Qty
13258	Stainless Steel Cylinder Stir Bars,	1000
	1.98mm x 4.80mm, for 2mL HPLC Vials	

#### **Magnetic Stir Bars**

Cat. No	. Description	Qty
50225	Magnetic Stir Bars, 5mm x 2mm, PTFE Coated	25
502100	Magnetic Stir Bars, 5mm x 2mm, PTFE Coated	100

#### **Accessories**

#### 12 x 32mm vials (sold separately)

Cat. No.	Description	Qty
11211-Case	11mm, 12 x 32, 2mL Wide Mouth Glass Crimp/Snap Vials	1000
11221-Case	11mm, 12 x 32, 2mL Wide Mouth Glass <b>Crimp</b> Vials	1000
11221M-Case	11mm, 12 x 32, 2mL Wide Mouth Glass Crimp Vials with Marking Spot	1000



## 48-Position Numbered Crimp Vial Assembly for Reaction Blocks

Cat. No.	Description	Qty
48221M	48-Position Numbered Crimp Vial	Each
	Assembly. Includes: Vials (11221M)	
	Marked 1-48; Stir Bar (13258) in Each	
	Vial; 48-Position Cap Mat (99948)	



Cap Mat for HPLC (2mL) Vials in 48-Position Block

Cat. No.	Description	Qty
99948	Clear Silicone/PTFE Cap Mat	5
	for HPC Vials in 48-Position	
	Reaction Block	









#### 1 DRAM, 24-Position Photoredox Catalysis Aluminum Reaction Blocks with 18mm Spacing, Gen II

1 DRAM (4mL, 15x45mm VIALS), GEN II

for 4mL (1 Dram), 15x45mm Vials

Catalog No.	Description	Qty
101240	Gen II, 24 Position, 5-BOLT, Photoredox Reaction Block for 15mm OD	Each
	(1 Dram) Vials, with 18mm Well Spacing	

## 1 DRAM, 24-Position Parallel Synthesis Aluminum Reaction Blocks with 18mm Spacing, Gen II for 4mL (1 Dram), 15x45mm Vials

Catalog No.	Description	Qty
104240	Gen II, 24 Position, 5-BOLT, Parallel Synthesis Reaction Block for 15mm	Each
	OD (1 Dram) Vials, with 18mm Well Spacing	







Replacement Parts (for photoredox AND parallel synthesis Gen II 24-position blocks)

Catalog No.	Description	Qty
101247	Top Rubber Mats for Gen II 24-Position Reactors (with 18mm Spacing)	25
101248	Top PFA Films for Gen II 24-Position Reactors (with 18mm Spacing)	25
SHS0114	Socket Head Screws for Gen II 24-Position Reactors and 96-Position Lightweight Reactors	5



Our 24-position Photoredox Catalysis Reactors with 18mm spacing (Gen II and original) are built specifically for use with our Lumidox® II 24-position, 18mm spaced Lens Mat Arrays

See page 24 for more information



analytical

## **Temperature Transfer Cover** for 24-Position, 18mm Spacing Reactors for 15mm OD Vials


Perfect for those who prefer to cap their vials instead of sealing them with the standard aluminum lid, rubber mat and PFA film that come with the block assembly. When using the block assembly in conjunction with a Thermal Transfer Deck+ (TTD+) and external liquid chiller, the cover helps control excessive heat generated by a Lumidox® II LED array and also greatly reduces light bleed.

The vials in the reactor block provide the alignment for the cover, but for extra stability users can secure the cover with four (4) bolts. The cover can also be flipped over and used as a reactor in itself or vial tray.

The Temperature Transfer Cover is available in a complete block assembly (101250) or as a separate part (249913).

Cat. No.	Description	Qty
101250	Gen II, 24-Position Photoredox Reaction Block (18mm Well Spacing) for 15x45mm (1 Dram) Vials, with Temperature Transfer Cover. Does NOT include rubber mat, PFA film and standard lid.	Ea
249913	Temperature Transfer Cover ONLY - for Gen II 24-Position Reactors with 18mm Well Spacing	Ea



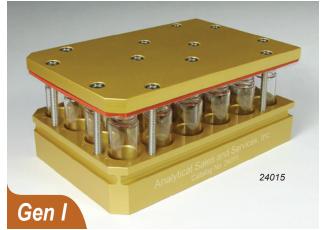

## Gen I



Use with 15 x 45mm vials, sold separately. See part number 31531-Case.



#### 31531

#### 1 DRAM, 24-Position Photoredox Catalysis Aluminum Reaction Blocks (20mm spacing)


for 4mL (1 Dram), 15x45mm Vials

Cat. No.	Description	Qty
24615	24-Position, Open Top/Open Bottom Aluminum Reaction Block for	Each
	15mm OD (1 Dram) Vials, Includes: Base Plate, Cover, Screws,	
	PFA Film and 2 Rubber Mats - for Mettler Toledo QX96	

## 1 DRAM, 24-Position Parallel Synthesis Aluminum Reaction Blocks (20mm spacing)

for 4mL (1 Dram), 15x45mm Vials

Cat. No.	Description	Qty
24015	24-Position Aluminum Reaction Block for 15mm OD (1 Dram) Vials, Includes: Base Plate, Cover, Screws, PFA Film and 2 Rubber Mats - for Mettler Toledo QX96	Each

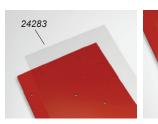






Replacement Parts (for photoredox AND parallel synthesis Gen II 24-position blocks)

Cat. No.	Description	Qty
24282	High-Temp Silicone/Rubber Cover for 24-Position Reaction Blocks for Vials	25
24283	PFA Film (0.005" Thick) for 24-Position Reaction Blocks for Vials	25
VSCREW24	1 3/4" Screws for 24-Position Aluminum Reaction Plate	100

#### 1 DRAM, 24-Position Photoredox Catalysis Aluminum Reaction Blocks with 18mm Spacing for 4mL (1 Dram), 15x45mm Vials

Cat. No.	Description	Qty
24626	24-Position Aluminum Reaction Block with 18mm spacing between well centers. For 15mm OD (1 Dram) Vials. Includes: Base Plate, Cover, Screws, PFA Film and 2 Rubber Mats	Each
24121	PFA Films for 24 Position, 18mm Spacing Reaction Blocks	25
24120	Silicone/Rubber Mats for 24 Position, 18mm Spacing Reaction Blocks	25
VSCREW24	1 3/4" Screws for 24-Position Aluminum Reaction Plate	100



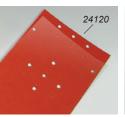



#### Vials and Caps

Cat. No.	Description	Qty
31531-Case	Advantage™ 13mm, 15 x 45, 4mL Clear Glass Screw Vials, Case	1000
31554-Case	PTFE Lined Cap	1000





## Photoredox Catalysis & Parallel Synthesis





## 2 DRAM, 24-Position Photoredox Catalysis Aluminum Reaction Blocks with 18mm Spacing, Gen II

For 17x60mm, 8mL (2 Dram) Glass Vials

Catalog No.	Description	Qty
103240	Gen II, 24 Position, 5-BOLT, Photoredox Reaction Block for 17mm OD	Each
	(2 Dram) Vials, with 18mm Well Spacing	

## 2 DRAM, 24-Position Parallel Synthesis Aluminum Reaction Blocks with 18mm Spacing, Gen II

For 17x60mm, 8mL (2 Dram) Glass Vials

Catalog No.	Description	Qty
106240	Gen II, 24 Position, 5-BOLT, Parallel Synthesis Reaction Block for 17mm OD (2 Dram) Vials, with 18mm Well Spacing	Each







Replacement Parts (for photoredox AND parallel synthesis Gen II 24-position blocks)

Catalog No.	Description	Qty
101247	Top Rubber Mats for Gen II 24-Position Reactors (with 18mm Spacing)	25
101248	Top PFA Films for Gen II 24-Position Reactors (with 18mm Spacing)	25
SHS6140	M6x1, 40mm Long, Low-Profile Socket Head Screw	5PK

#### Gen II / Gen II

#### **NEW!** Para-dox® Gen II Aluminum Reaction Blocks

#### For Photoredox Catalysis and Parallel Synthesis

#### Benefits of the Gen II line include:

- Less bolts = less time to assemble
- Less light bleed when used with Lumidox<sup>®</sup> II LED arrays (48 and 96 position blocks)
- Less Maintenance
- Vials fit tighter (less shifting)
- Longer lasting screws (larger diameter, larger threads)

#### **Additional features:**

- Port for thermocouple, sensors, etc.
- Corner holes on both the lid and the bottom for accessory attachment
- · Larger holes in lid to accommodate a variety of needle sizes
- Uses only one 1/8" rubber mat compared to two 1/16" mats on previous versions (makes for easier assembly)

### Parallel Synthesis vs. Photoredox Catalysis Reaction Blocks

Reactors for both applications are nearly identical. The difference lies in the base (vial rack) of the block. Reactors for parallel synthesis have wells with closed bottoms, whereas photoredox blocks have wells with open holes in the bottom allowing for light transmission via a Lumidox®II Array, as well as to read 2D barcodes.





Photoredox Catalysis Block, Gen II









Use with 17 x 60mm vials, sold separately. See part number 31760-Case.



31760



## 2 DRAM, 24-Position Photoredox Catalysis Aluminum Reaction Blocks


for 8mL (2 Dram), 17x60mm Vials

Cat. No.	Description	Qty
24617	24-Position, <b>Open Top/Open Bottom</b> Aluminum Reaction Block for 17mm OD (2 Dram) Vials, Includes: Base Plate, Cover, Screws, PFA Film	Each
	and 2 Rubber Mats	

## 2 DRAM, 24-Position Parallel Synthesis Aluminum Reaction Blocks

for 8mL (2 Dram), 17x60mm Vials

Cat. No.	Description	Qty
24017	24-Position Aluminum Reaction Block for 17mm OD (2 Dram) Vials.	Each
	Includes: Base Plate, Cover, Screws, PFA Film and 2 Rubber Mats	

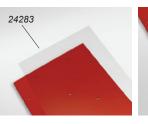





Replacement Parts (for photoredox AND parallel synthesis Gen II 24-position blocks)

Cat. No.	Description	Qty
24282	High-Temp Silicone/Rubber Cover for 24-Position Reaction Blocks for Vials	25
24283	PFA Film (0.005" Thick) for 24-Position Reaction Blocks for Vials	25
VSCREW24	1 3/4" Screws for 24-Position Aluminum Reaction Plate	100

#### Vials and Caps

Cat. No.	Description	Qty
31760-Case	Advantage™ 17 x 60mm, 8mL Clear Glass Screw Vials, Case	1000
31542-CASE	15mm Solid Black Polypropylene Screw Caps with PTFE/F217 Liners	1000
31543-CASE	Black Open Top PP Cap with PTFE/Silicone (0.065") Setpa, 15-425mm Thread	1000





#### Accessories

Cat. No.	Description	Qty
84001-Case	1mL Clear Glass Shell Vials, 8 x 30mm	1000
13258	Stainless Steel Cylinder Stir Bars, 1.98mm x 4.80mm (for 8x30mm Vials)	1000
SD1000	Milwaukee M4 1/4in. Hex Screwdriver Kit. Includes 2 Batteries and Charger (Note: batteries not available for international shipping)	Each













#### **Lumidox® II Line of products:**

- 96-position LED Arrays
- 48-position LED Arrays for TCR (see page 4)
- 24-position Arrays
- LumLamps
- Lumidox II Controller
- Cell Culture Plate Alignment Spacers for Lumidox II LED Arrays (see website)

#### **Key Features:**

- Reduces photochemical reaction time
- Up to 5 selectable output STAGES
- Up to 15 wavelengths available
- Higher optical/radiometric power than previous Lumidox line of products
- Patented LED array design
- Experiment with electro-photochemical reactions by pairing with HTe<sup>-</sup>Chem Electrochemistry Assemblies (see page 30)



## For Photoredox Catalysis Applications

Taking advantage of recent breakthroughs in LED technology, Analytical now offers the next generation Lumidox® II with higher optical/radiometric power than ever before. Lumidox II offers a multitude of unique and complimenting photonic devices, in varying wavelengths, amplitudes and footprints.

The Lumidox II collection includes new LED Arrays with a wider range of options to choose. Also, check out our LumLamp for experimental applications. All of our Lumidox II illumination devices are controlled by the Lumidox II controller.

## Lumidox II Arrays and LumLamps are available in up to 15 wavelengths:

365 (UV365) 405 (UV405) 505 (Cyan) 660 (Deep Red) 375 (UV375) 420 (Violet) 527 (Green) 730-IR (Infrared) 385 (UV385) 445 (Indigo) 590 (Amber) White 395 (UV395) 470 (Blue) 630 (Red)

Analytical has also developed a myriad of complimentary apparatus that works with Lumidox. These range from SLAS cooling blocks to our Temperature Controlled Reactors (TCR, pages 4, 5). We also offer high throughput electrochemical assemblies that can work with Lumidox to expand your experimentation into electro-photochemical reactions (see page 30).

# Lumidox II Controller Lumidox II PUSH ON INTERPRETATION FORCE 9, 1 H ENDERNOT FORCE 9, 1 H

#### **Peripheral Devices**





#### Lumidox® II Controller

- Controls the output of a Lumidox Array or LumLamp
- Select one of 5 levels of optical power (STAGES).
   STAGES are displayed in easy to read, easy to calculate whole numbers.
- Onboard illumination elapsed timer
- Automatic countdown shut-off timer
- USB Enabled API for advanced users running their setups remotely

Catalog No.	Description
LUM2CON	Lumidox II LED Controller. Includes: Controller,
	Power Supply, Power Cable, Manual



#### PRODUCT NOTE:

Arrays and LumLamps are tuned in-house. Calibration data is stored onboard (not on the controller), allowing the use of different illuminators with one controller. Note: the controller can operate different Lumidox devices, but only 1 device at a time.

Controller accepts 100-240 voltage – refer to the Lumidox® II manual for information regarding usage, electrical requirements, etc.

#### **LED Output STAGES**

By default, all Lumidox II devices are factory calibrated with 5 discrete linearly stepped output STAGES. STAGE 1 output is the least radiometric power while STAGE 5 output is the most. STAGES are calibrated to the nearest whole number of radiant flux, and displayed in milliwatts (mW).

Linear Step (default)

STAGE 1 - 50mW STAGE 2 - 100mW STAGE 3 - 150mW

STAGE 4 - 200mW STAGE 5 - 250mW



#### **Custom LED Tuned STAGES**

Analytical offers custom tuned stages to meet specific and unique requirements. These can range from maximum light output to low output to tight resolution.

**NOTE:** STAGE settings are stored on each individual Lumidox® II **peripheral** device (LED Array or Lumlamp), and are **not tunable via the controller**. Custom STAGE settings can only be calibrated by Analytical Sales and Services prior to shipment. Please contact Analytical before ordering.

**High Output** (for more aggressive reactions)

STAGE 1 - 100mW STAGE 2 - 200mW STAGE 3 - 400mW

STAGE 4 - 600mW STAGE 5 - 800mW High Resolution

(for greater precision)

STAGE 1 - 100mW


STAGE 2 - 110mW STAGE 3 - 115mW

STAGE 4 - 120mW STAGE 5 - 125mW **Low Output** (good for cell culture and PCR)

STAGE 1 - 10mW STAGE 2 - 15mW STAGE 3 - 20mW STAGE 4 - 25mW STAGE 5 - 50mW





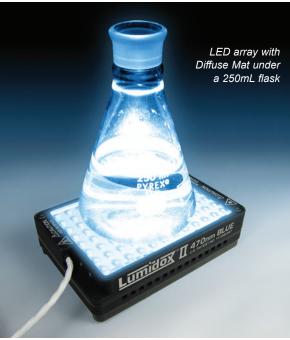

#### **Lumidox® II LED Arrays**

U.S. Patent No. 11,458,447

Our patented LED arrays provide wavelength and power specific illumination to samples for photoredox catalysis applications.

- Up to 15 different color wavelengths are available\*
- 5 output STAGES (customizeable) with calibration data stored on board\*\*
- Available in two Surface Mat styles
- Three options for Base configurations

#### Radiant Flux Values per STAGE - Example

(typical values for 96-Position LED Arrays with Active Cooling Base and Lens Mat)

	STAG	E 1	STAG	E 2	STAG	E 3	STAG	E 4	STAG	E 5
Wavelength	Per well (mW)	Total (W)								
UV365	25	2.4	55	5.3	80	7.7	105	10.1	135	13.0
UV375	25	2.4	50	4.8	75	7.2	110	10.6	140	13.4
UV385	30	2.9	65	6.2	100	9.6	145	13.9	165	15.8
UV395	30	2.9	65	6.2	105	10.1	140	13.4	170	16.3
UV405	25	2.4	60	5.8	90	8.6	125	12	160	15.4
420-Violet	30	2.9	60	5.8	90	8.6	120	11.5	150	14.4
445-Indigo	60	5.8	120	11.5	195	18.7	240	23.0	295	28.3
470-Blue	45	4.3	95	9.1	140	13.4	180	17.3	220	21.1
505-Cyan	35	3.4	65	6.2	90	8.6	115	11.0	135	13.0
527-Green	25	2.4	55	5.3	80	7.7	100	9.6	110	10.6
590-Amber	75	7.2	95	9.1	115	11.0	150	14.4	180	17.3
630-Red	30	2.9	55	5.3	85	8.2	115	11.0	145	13.9
660-Deep Red	40	3.8	70	6.7	105	10.1	140	13.4	170	16.3
IR730	35	3.4	65	6.2	95	9.1	125	12.0	155	14.9
White	50	4.8	100	9.6	150	14.4	200	19.2	300	28.8

Irradiance charts can be found on our website

#### **Surface Mat Styles**

Array top surfaces come in two styles, Lens Mat and Diffuse Mat.



#### **Lens Mat Surface**

- Ultra-clear, molded to fit into the holes of a Para-dox® Reaction Block
- Captures nearly all light emitted by the array's LEDs and directs it into the vials in the reaction block
- · Chemically inert silicone



#### **Diffuse Mat Surface**

- Flat surface, making for easy pairing with SLAS footprint apparatus
- Can be used for niche applications such as illuminating cell culture flasks, reservoir plates, large scale containers, etc.
- Chemically inert silicone


#### **About Array Cooling and Base Options**

Lumidox® II LED Arrays can generate a considerable amount of heat at any output stage and therefore need to be cooled. Both **Lens Mat and Diffuse Mat** style arrays are available with an **Active Cooling** base, **Solid** base or **Flow-Through** base. Each has it's own specific way of cooling the LEDs in the array and can affect different application situations. *Note: solid and flow-through base arrays are not self-cooling and require the use of a sufficient external chiller/cooling device.* 

#### **Array Base Comparison**

#### **Active Cooling Base**



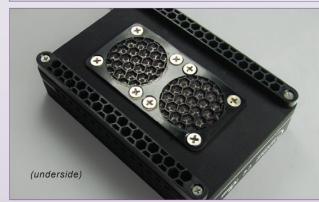

- Self-cooling, no external cooling source required
- Fully conforms to SLAS/ANSI standard dimensions
- Ideal for lower output applications like cell culture and PCR work (may require special Alignment Spacer)
- Not compatible with tumble stirrers or devices that generate powerful magnetic fields
- · Can be placed in an incubator
- \* For comparison, Radiant Flux Value charts for Active Cooling Base and Solid/Flow-through Base Arrays are posted on our website.

Note: All LED arrays can operate in an incubator at 37°C, 95% humidity

#### **Solid Base**





External cooling source is required (a Thermal Transfer Deck with a recirculating liquid chiller is recommended)

- A Thermal Transfer Deck (TTD, see page 25) can be attached to the base when used in conjunction with a recirculating chiller. The TTD can be removed if using a different cooling source, such as a cooling bay
- Fully conforms to SLAS/ANSI standards and can be used with cooling baths, plates, or other cooling chamber (if NOT being used with our Thermal Transfer Deck and recirculating liquid chiller)
- Ideal for high output requirements up to nearly 3x more radiometric power output than Active Base arrays\*
- · Best highest output, highest intensities



#### Flow-Through Base

Recirculating chiller required, liquid flows through the LED array base

- Not recommended to be used with the Temperature Controlled Reactors
- Nearly 3x the radiometric power output of Active Cooling Base arrays
- Short & lightweight improved usability with orbital shakers due to lower overall mass



<sup>\*</sup> Most LED arrays (excluding Discovery, page 23) are wavelength-specific - only one wavelength per device

<sup>\*\*</sup> Calibration data is stored onboard (not on the controller), allowing the use of different illuminators with one controller. Controller can only operate 1 device at a time. STAGE settings are NOT user-adjustable and must be configured prior to shipping





#### 96-Position LED Arrays







Solid Base Flow-Through Base

96-Position LED Arrays with <b>Lens Mat</b>		<b>Active Cooling Base</b>	Flow-Through Base*	Solid Base*
Wavelength (nm)	Description	Catalog No.	Catalog No.	Catalog No.
UV365	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA365	LUM296LF365	LUM296LS365
UV375	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA375	LUM296LF375	LUM296LS375
UV385	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA385	LUM296LF385	LUM296LS385
UV395	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA395	LUM296LF395	LUM296LS395
UV405	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA405	LUM296LF405	LUM296LS405
420-VIOLET	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA420	LUM296LF420	LUM296LS420
445-INDIGO	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA445	LUM296LF445	LUM296LS445
470-BLUE	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA470	LUM296LF470	LUM296LS470
505-CYAN	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA505	LUM296LF505	LUM296LS505
527-GREEN	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA527	LUM296LF527	LUM296LS527
590-AMBER	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA590	LUM296LF590	LUM296LS590
630-RED	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA630	LUM296LF630	LUM296LS630
660-DEEP RED	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA660	LUM296LF660	LUM296LS660
730-IR	Lumidox II 96-Position LED Array with Lens Mat	LUM296LA730	LUM296LF730	LUM296LS730
WHITE	Lumidox II 96-Position LED Array with Lens Mat	LUM296LAWHT	LUM296LFWHT	LUM296LSWHT

\* requires the use of a sufficient chiller/cooling device

See pages 20-21 for details on surface mats and bases









Flow-Through Base

96-Position LED Arrays with <b>Diffuse Mat</b>		<b>Active Cooling Base</b>	Flow-Through Base*	Solid Base*	
Wavelength (nm)	Description	Catalog No.	Catalog No.	Catalog No.	
UV365	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA365	LUM296DF365	LUM296DS365	
JV375	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA375	LUM296DF375	LUM296DS375	
JV385	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA385	LUM296DF385	LUM296DS385	
UV395	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA395	LUM296DF395	LUM296DS395	
JV405	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA405	LUM296DF405	LUM296DS405	
120-VIOLET	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA420	LUM296DF420	LUM296DS420	
145-INDIGO	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA445	LUM296DF445	LUM296DS445	
170-BLUE	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA470	LUM296DF470	LUM296DS470	
505-CYAN	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA505	LUM296DF505	LUM296DS505	
527-GREEN	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA527	LUM296DF527	LUM296DS527	
590-AMBER	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA590	LUM296DF590	LUM296DS590	
30-RED	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA630	LUM296DF630	LUM296DS630	
660-DEEP RED	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA660	LUM296DF660	LUM296DS660	
730-IR	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DA730	LUM296DF730	LUM296DS730	
VHITE	Lumidox II 96-Position LED Array with Diffuse Mat	LUM296DAWHT	LUM296DFWHT	LUM296DSWHT	



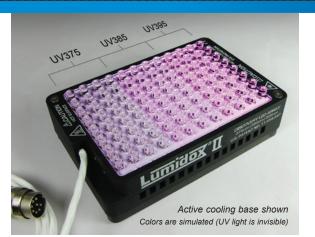

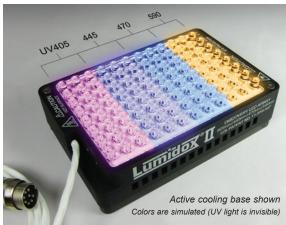





#### **Discovery/Screening 96-Position LED Arrays**

Lumidox® II Discovery LED Arrays for screening offer scientists an economical and efficient way to evaluate most of the wavelengths that the Lumidox II line of products has to offer. Discovery LED arrays come in varying wavelength combinations and are available with an active cooling, solid or flow-through base.

There are currently 15 wavelength options available in the Lumidox II product line. With Discovery LED arrays, 9 of those wavelengths are combined into 3 units (no need to purchase 9 individual units). The three models of Discovery LED arrays come in these wavelength combinations:

- Discovery 1 (3 UVs):
   375nm, 385nm, 395nm (4 columns of 8 LEDs / wavelength)
- Discovery 2 (UV405, Indigo, Blue & Amber):
   405nm, 445nm, 470nm, 590nm (3 columns of 8 LEDs / wavelength)
- Discovery 3 (Cyan, Green): 505nm, 527nm (6 columns of 8 LEDs / wavelength)

The six remaining wavelengths in the Lumidox II line - UV365, 420-VIOLET, 630-RED, 660-DEEP RED, 730-IR, and White - are not compatible for placement in Discovery LED arrays due to varying electrical conditions. They can, however, be purchased as single wavelength units to complete your LED array collection.

#### 96-Position Discovery LED Arrays with Lens Mat

		Active Cooling Base	Flow-Through Base*	Solid Base*
Wavelength (nm)	Description	Catalog No.	Catalog No.	Catalog No.
UV375 UV385 UV395	Discovery 1 LED Array w/ Lens Mat (4 rows of 8 LEDs per wavelength)	LUM296LAG1	LUM296LFG1	LUM296LSG1
UV405				
445-INDIGO 470-BLUE	Discovery 2 LED Array w/ Lens Mat (3 rows of 8 LEDs per wavelength)	LUM296LAG2	LUM296LFG2	LUM296LSG2
590-AMBER				
505-CYAN 527-GREEN	Discovery 3 LED Array w/ Lens Mat (6 rows of 8 LEDs per wavelength)	LUM296LAG3	LUM296LFG3	LUM296LSG3
Single Wavelengt	th LED Arrays			
UV365	96-Position LED Array w/ Lens Mat	LUM296LA365	LUM296LF365	LUM296LS365
420-VIOLET	96-Position LED Array w/ Lens Mat	LUM296LA420	LUM296LF420	LUM296LS420
630-RED	96-Position LED Array w/ Lens Mat	LUM296LA630	LUM296LF630	LUM296LS630
660-DEEP RED	96-Position LED Array w/ Lens Mat	LUM296LA660	LUM296LF660	LUM296LS660
730-IR	96-Position LED Array w/ Lens Mat	LUM296LA730	LUM296LF730	LUM296LS730
WHITE	96-Position LED Array w/ Lens Mat	LUM296LAWHT	LUM296LFWHT	LUM296LSWH
	·			

<sup>\*</sup> requires the use of a sufficient chiller/cooling device

#### **Thermal Transfer Decks**

See page 25 for more information

Cat. No.	Description
266530	SLAS Thermal Transfer Deck (TTD)
266540	TTD+ (Thermal Transfer Deck with High-end Fittings)
266550	SLAS Footprint, "Straight-Thru" Thermal Transfer Deck













#### Lumidox II 24-Position LED Arrays are available in the these wavelengths:

365 (UV365)	395 (UV395)	470 (Blue)	590 (Amber
375 (UV375)	405 (UV405)	505 (Cyan)	630 (Red)
385 (UV385)	445 (Indigo)	527 (Green)	White

#### Lumidox® II 24-position LED Arrays with 9mm Spacing

- Typically used with a Para-dox® 24-well Reaction Block for 8x30mm vial inserts (see pages 10, 11)
- Can also be used with a HTe-Chem Electrochemistry Assembly (see page 30)

		<b>Active Cooling Base</b>	Solid Base*
Wavelength (nm)	Description	Catalog No.	Catalog No.
UV365	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA365	LUM224LS365
JV375	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA375	LUM224LS375
JV385	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA385	LUM224LS385
JV395	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA395	LUM224LS395
JV405	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA405	LUM224LS405
145-INDIGO	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA445	LUM224LS445
170-BLUE	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA470	LUM224LS470
505-CYAN	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA505	LUM224LS505
527-GREEN	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA527	LUM224LS527
590-AMBER	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA590	LUM224LS590
30-RED	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LA630	LUM224LS630
WHITE	24-Well LED Array with 9mm Spacing, Lens Mat	LUM224LAWHT	LUM224LSWHT

\* requires the use of a sufficient chiller/cooling device

#### See pages 20-21 for details on surface mats and bases





#### **Lumidox® II 24-position LED Arrays with 18mm Spacing**

Built specifically for use with our Para-dox® 24-well, 18mm spaced **Reactor Blocks** 

See pages 14, 16 for more information

		<b>Active Cooling Base</b>	Solid Base*
Wavelength (nm)	Description	Catalog No.	Catalog No.
UVC-255	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA255	LUM22418LS255
UV365	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA365	LUM22418LS365
UV375	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA375	LUM22418LS375
UV385	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA385	LUM22418LS385
UV395	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA395	LUM22418LS395
UV405	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA405	LUM22418LS405
420-VIOLET	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA420	LUM22418LS420
445-INDIGO	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA445	LUM22418LS445
470-BLUE	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA470	LUM22418LS470
505-CYAN	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA505	LUM22418LS505
527-GREEN	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA527	LUM22418LS527
590-AMBER	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA590	LUM22418LS590
630-RED	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LA630	LUM22418LS630
WHITE	24-Well LED Array with 18mm Spacing, Lens Mat	LUM22418LAWHT	LUM22418LSWHT

analytical





#### TECH NOTE:

If using in conjunction with a Flow Reactor (page 28), a Thermal Transfer Deck with High-end Fittings (TTD+) must be used



Cat. No.	Description	Qty
266540	TTD+ (Thermal Transfer Deck with High-end Fittings)	Each
266550	SLAS Footprint, "Straight-Thru" Thermal Transfer Deck+	Each
	(Thermal Transfer Deck with High-end Fittings)	

#### **Thermal Transfer Decks**

- · Must be used with a recirculating liquid chiller or heater
- Can be attached to a Lumidox® II solid base LED arrays for direct cooling of the array

When connected to an external recirculating liquid chiller or heater, our Thermal Transfer Decks can be used as both chilling plates or hot plates. Apply heat or cold directly to your samples or sample vessels (such as Para-dox® Reaction Blocks), or use it to cool your Lumidox® II solid base LED array.

Thermal Transfer Decks may be linked in novel ways to form gradients, or to serve other more experimental needs. Push-toconnect fittings allow for quick tubing insertion/removal. Input and Output ports are user configurable, and may be relocated between side and top faces\*. Units are shipped with side exit ports as standard. Decks are stackable, and multiple decks can be chained together for a total temperature control solution. Compatible with a wide variety of coolants (water, propylene glycol, ethylene glycol, etc.).

\* Ports are not user configurable on the Straight-Thru TTD




Attach a Thermal Transfer Deck+ to a Lumidox II SOLID BASE LED Array for direct cooling of the array



Use an additional TTD+ to transfer heat or cold directly to your reaction block samples





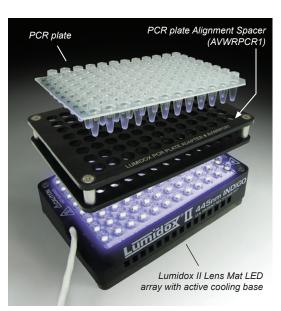

**Lumidox**®**II** 

Lumidox / PCR Setup



Lumidox / Cell Culture Setup



#### Lumidox® II for Cell Culture / Biological Applications

We can support your photon irradiation needs for your cell culture or biological applications, such as:

- Photoconverting live mammalian cells with Dendra protein for fluorescent imaging days after photoconversion
- Testing of light-induced toxicity of certain compounds on different cell lines for photodynamic therapy applications
- Studying the photothermal effects of compounds in live cells
- Testing photodynamic therapy of various innovative compounds on live cells
- The use of photodynamic therapy as a type of cancer treatment

#### **Lumidox II LED Arrays with Active Cooling Base**

Lumidox II LED arrays generate wavelength specific light for your cell culture or biological experimentation.

- Available in 15 wavelengths from UV, Visible through IR.
- Run at 37°C indefinitely (70°C max) at humidity <=99%.
- Active Cooling Base LED arrays are fully compatible with incubation chambers.
- Use Lens Mat Arrays for 24 or 96 well flat bottom collection plates, or with Analytical's 48-well TCR.
- Use **Diffused Mat** Arrays for cell culture flasks, petri dishes, larger 12-well and 6-well cell culture plates, etc..
- Active Cooling Base arrays are recommended for bio/cell culture work, but other base configurations are available.

#### **Cell Culture Plate Alignment Spacers**

For Low Output (<50mW) Experimentation

Designed for low output applications using **Lens Mat** arrays (not compatible with diffuse mat arrays). Run your array at a stage with an appropriate output and environment. Higher output levels and temperatures can deform the alignment spacer. While these Alignment Spacers provide the most practical method for consistent positioning of your samples over the LEDs, if higher output (>50mW per LED) is desired, you can manually position your sample plate over the array without a spacer.

- All our Low Profile collection plates
- Cellvis P96-1.5H-N
- Greiner 655180
- · Corning Costar cell culture plates
- Eppendorf cell culture plates (contact us with your specific plate)
- Thermo Scientific Nunc MicroWell plates
- Perkin Elmer ViewPlate (formerly Packard ViewPlate)

Alignment Spacers are available for any standard cell culture plate by request.

#### Lumidox® II Controller

- Five stages of light intensity to choose from that can be calibrated to your custom levels at the time of purchase see STAGE settings, page 19.
- Controller has an onboard timer to reproducibly limit exposure.









**24 well setup** with Lumidox 24-well LED Array - See page 24 for array options

#### Lumidox® II LED Arrays with Active Cooling Base

Recommended wavelengths for biological and cell culture work. Other wavelengths are available.

Catalog No.	Description	Wavelength
LUM296LA445	Lumidox® II 96-position LED Array with Lens Mat & Active Cooling Base	445nm - Indigo
LUM296LA470	Lumidox® II 96-position LED Array with Lens Mat & Active Cooling Base	470nm - Blue
LUM296LA527	Lumidox® II 96-position LED Array with Lens Mat & Active Cooling Base	527nm - Green
LUM296LA630	Lumidox® II 96-position LED Array with Lens Mat & Active Cooling Base	630nm - Red
LUM296LA660	Lumidox® II 96-position LED Array with Lens Mat & Active Cooling Base	660nm - Deep Red
LUM296LAWHT	Lumidox® II 96-position LED Array with Lens Mat & Active Cooling Base	White

#### Cell Culture Plate Alignment Spacer

ocii oait	on Sultare Flate Anginient Spacer		
Catalog No.	Description		
A655180	Alignment Spacer for Lumidox II LED Arrays, for use with Greiner 655180 Plate		
AP9615HN	Alignment Spacer for Lumidox II LED Arrays, for use with Cellvis P96-15H-N Well Plates		
A204625	Alignment Spacer for Lumidox II LED Arrays, for use with Cell Culture Plate Part # 96624		
A204627	Alignment Spacer for Lumidox II LED Arrays, for use with Cell Culture Plate Part # 96626		
ACRCLP01	Alignment Spacer for Lumidox II LED Arrays, for use with Corning Cell Culture Plates		
A6005430	Alignment Spacer for Lumidox II LED Arrays, for use with Perkin Elmer ViewPlate (Packard ViewPlate)		
A204628	Alignment Spacer for Lumidox II LED Arrays, for use with Cell Culture Plate Part # 384628		
ACRCLP03	Alignment Spacer for Lumidox II LED Arrays, for use with Corning 2592 Plate		
ACRCLP02	Alignment Spacer for Lumidox II LED Arrays, for use with Corning 3631 Plate		
A655097	Alignment Spacer for Lumidox II LED Arrays, for use with Greiner 655097 Plate		
AT11028	Alignment Spacer for Lumidox II 24-position LED Arrays, for use with 24 Deep Well Plate #24128		

#### **Cell Culture Plates**

Catalog No.	Description	QTY
96624	96-well Cell Culture/Imaging Microplate (Agilent) - Clear with Flat Bottom, 400µL Per Well	40
96518	96 Well, Flat Bottom Clear PS Microplate, 25-340μL working volume per well	100
96626	96-well Cell Culture/Imaging Microplate (Agilent) - Black with Clear Flat Bottom, 400µL Per Well	32
961515	Cellvis 96-well Black Polystyrene/Glass Bottom Plate with high performance #1.5 cover glass	20
96097	$96 \ Well, \ Black/Clear \ Flat \ Bottom \ Well \ PS \ Microplate, \\ 25-340 \mu L \ working \ volume \ per \ well, \ high \ binding, \ sterile$	40
384628	384-well Cell Culture/Imaging Microplate, 140µL Per Well	32
24128	24 Deep Well Plate, V-Bottom	50

Please see our website for additional collection plates



#### **PCR Plate Alignment Spacer**


Catalog No.	Description
AVWRPCR1	PCR Plate Alignment Spacer for Lumidox®II 96-position LED Lens Mat Arrays


#### **Borofloat 33 Glass Block**

Use to reduce the amount of light intensity delivered to your samples by increasing the distance between the samples and the light source (LED array).

`	• /
Catalog No.	Description
SLASBG33	Borofloat 33 Glass Block



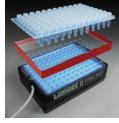

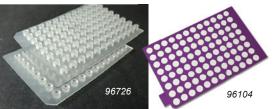

#### Cap Mats and Films for PCR Plates

Cat. No.	Description	QTY
96726	6mm Silicone/PTFE Cap Mats for 96-well PCR Collection Plates	5
96104	Pattern Adhesive <i>Resealable</i> Silicone Sealing Film, Round 96-Well Pattern	5







**Lumidox**® **II** 

#### Flow Reactor

The Para-dox® Flow Reactor offers an alternative way to conduct photoreactive research on fluidic samples. Used in conjunction with a Lumidox®II Diffuse Mat Array, wavelength-specific light can be applied to sample fluid circulating through 1/8" OD tubing. The Flow Reactor can be liquid cooled or heated and is highly configurable. Add a Thermal Transfer Deck for a complete flow solution.

Users may configure the Flow Reactor in various ways, including multiple input streams using the machined channels. Tubing can be replaced and installed by the user with a tubing tamp tool.

- SLAS/SBS footprint
- Dripless CPC fittings
- 1/8" OD unions for ease of analyte injection
- User-configurable output/input locations, side or top (factory set to side).
- Exposure volume of 2mL using 1/8" OD, 1/16" ID tubing.
   The exposure volume may be set to 1mL if two streams are used.
- FEP Tubing Provided
- Cradled design reduces light bleed and helps "seat" and align Lumidox unit
- Compatible with water, glycol based coolants, and silicone based coolants

#### Flow Reactor

Catalog No.	Description	Qty
266100	Flow Reactor	Each





#### Recommended Chillers:

Cat. No.	Description
TCube Edge	TCube Edge Recirculating Chiller
MC280OLE	Minichiller 280 Ole Recirculating Chiller













## A lamp system specifically designed for your experimental needs!

#### LumLamp

The LumLamp system stands apart with multiple wavelengths, tight linewidths, and remote (corded) operation. It has radiant flux levels up to 3W. LumLamps are directly controlled with our Lumidox® II Controller, providing optical stability, user configurable timing and output configuration.

With 4 differing reflector types, 3 styles of lens, and up to 12 wavelengths, LumLamp is available in 104+\* possible configurations.

#### Lens Types:







#### **Reflector Types:**

15° Narrow Spot Beam 30° Medium Beam 40° Wide Beam 80° Extra Wide Beam

#### LumLamps are available in the following wavelengths:

365 (UV365)	395 (UV395)	470 (Blue)	590 (Amber
375 (UV375)	405 (UV405)	505 (Cyan)	630 (Red)
385 (UV385)	445 (Indigo)	527 (Green)	White

<sup>\*</sup> Contact us for custom wavelengths

#### Lumidox® II Controller

- Controls the output of a Lumidox Array or LumLamp
- 5 individually calibrated levels of optical power (STAGES)
- Optical power levels are provided in easy to read, easy to calculate whole numbers
- Onboard illumination elapsed timer
- Automatic countdown shut off timer
- USB Enabled API for advanced users running their setups remotely

#### **Lumidox II LED Controller**

Catalog No.	Description	Qty
LUM2CON	Lumidox II LED Controller.	Each
	Includes: Controller, Power Supply, Power Cable, Manual	



Lumidox Controllers come with one pair of UV protective goggles. Additional goggles can be purchased.

#### **UV Safety Goggles**


-		
Catalog No.	Description	Qty
UVGoggles	Additional UV Safety Goggles	Each





Electrochemistry HTe<sup>-</sup>Chem Assemblies Electrochemistry



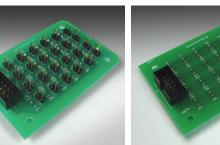



#### HTe-Chem Electrochemistry Assemblies

Analytical Sales' new **HTe<sup>-</sup>Chem** line of products allows for broad electrochemical reactions (such as electrosynthesis, organic electrochemistry, electrophotochemistry, etc.) to be carried out on Analytical Sales standard 24-well minirack platform. HTe<sup>-</sup>Chem allows you to accelerate your electrochemistry workflow by permitting multiple simultaneous constant current, constant voltage, and/or electrophotochemical experiments to be run.

There are two optimized setups to choose from:

The **constant current** setup allows the user to set a precise current, which will be maintained by the power supply throughout the reaction. The voltage will automatically be varied based on the changing electrical conditions as the reaction progresses.

The **constant voltage** setup allows the user to set a precise voltage, which will be maintained by the power supply throughout the reaction. In this setup, the current will automatically be varied based on the changing electrical conditions as the reaction progresses.

#### HTe<sup>-</sup>Chem Assemblies & Kits

Cat. No.	Description	Qty
700100	Constant Current Electrochemistry Assembly	Each
700150	Constant Current Electrochemistry Kit (Includes Electrode Kit)	Each
700200	Constant Voltage Electrochemistry Assembly	Each
700250	Constant Voltage Electrochemistry Kit (Includes Electrode Kit)	Each

#### Electrodes for HTe<sup>-</sup>Chem Assemblies

Analytical Sales offers a myriad of electrode choices which can be used as either cathodes or anodes depending on your experimental needs.


Cat. No.	Description	Qty
700500	Graphite Electrodes	30
700500-250	Graphite Electrodes - Bulk Pack	250
700525	Lead Electrodes	25
700550	Tin Electrodes	25
700575	Cadmium Electrodes	25
700600	Zinc Electrodes	25
700600-250	Zinc Electrodes - Bulk Pack	250
700650	Platinum Electrodes	25
700656	Platinum Electrodes - Row Pack	6
700675	Iron Electrodes	25
700700	Stainless Steel Electrodes	25
700700-250	Stainless Steel Electrodes - Bulk Pack	250
700725	Cobalt Electrodes	25
700750	Nickel Electrodes	25
700800	NiChrom Electrodes	25
700800-250	NiChrom Electrodes - Bulk Pack	250
700825	Silver Electrodes	25
700850	Copper Electrodes	25
700875	Tungsten Electrodes	25
700900	Aluminum Electrodes	25
700950	Magnesium Electrodes	25
700400	Electrode Kit. Includes One Pack of Each of Graphite, Zinc, Stainless Steel, Nickel, NiChrom, Aluminum, Copper and Magnesium Electrodes (platinum not included)	Each









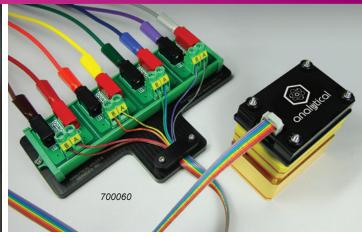












#### **Power Supply and Accessories**

Cat. No.	Description	Qty
700050	Calibrated DC Power Supply, 4 Output Multi Range, 420W Combined Output	Each
700060	Breakout Cable Assembly, for use with Calibrated DC Power Supply	Each
700070	Ribbon Cable (if using alternate power supply)	Each
701020	Replacement PCB Jumpers (for constant current setup)	100
701015	Upper Gasket (recommended: replace after 3-5 uses)	5
701005	Lower Gasket (recommended: replace after 3-5 uses)	5
701025	PFA Sheet for HTe <sup>-</sup> Chem Assembly (recommended: replace after every use)	25
24260	Bottom Rubber Mats	25
84001-CASE	1mL Clear Glass Shell Vial	1000

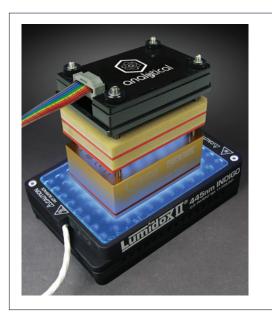



For experimentation with electrophotochemical reactions, HTe<sup>-</sup>Chem assemblies are fully compatible for use with Lumidox<sup>®</sup>II 24-well, 9mm spaced LED arrays.

See page 24 for more information




**Electrochemistry Electrochemistry Flow Electrolysis** Flow Electrolysis

#### Flow Electrolysis

Analytical is pleased to offer a solution for electrolysis experimentation in the form of a standardized flow electrolysis setup. This new and upcoming technology allows for rapid and efficient production of complex chemical compounds via flow electrolysis. With an interelectrode gap smaller than a millimeter, and the introduction of flow geometries that promote turbulent mixing, our Flow Electrolysis Cell allows for heightened selectivity and yield when compared to flask based reactors. Additionally, the Flow Electrolysis Cell requires minimal new equipment to use. With fittings for 1/8" OD HPLC tubing, fluid flow and electrodes can be connected to power supplies via standard connectors will automatically be varied based on the changing electrical conditions as the reaction progresses.

#### **Undivided Cell (single stream)**

In an undivided cell, both electrodes are washed with the same electrolyte and only one fluid circuit is used; the opposite side of the cell is sealed with HPLC plugs and an electrode without through-holes. This allows for low-resistance reactions which will output combined products.

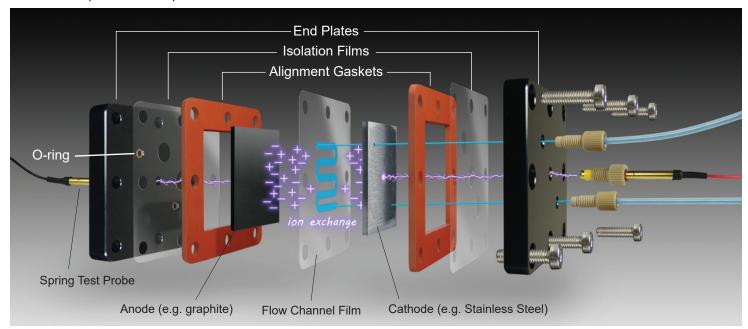





#### Flow Electrolysis Undivided Cell (single stream) Assembly Assembly includes:

- Impervious Graphite Plate Electrode (anode)
- 316L Stainless Steel Plate Electrode (cathode)
- Flow Channel PFA Films (1 Meandering, 1 Tangential, 1 Simple,
- 1 Fin Separator, 1 Flooded Flow)
- Isolation Layer PFA Films (2)
- Orange Silicone Rubber Alignment Gaskets (2)
- Viton<sup>™</sup> Chemical-resistant Internal O-rings, 2mm ID (2)
- Viton<sup>™</sup> Chemical-resistant External O-rings, 3mm ID (2)
- Spring Test Probes/Pogo Pins (1 red, 1 black)
- Aluminum End Plates (2)
- · Fittings and Hardware

at. No.	Description	Qty
1088100	Flow Electrolysis Undivided Cell Assembly (single stream)	Each

#### **Channel Film Volume**

Film Type	Surface Area (mm^2)	Volume at 0.254mm Thick (mm^3)	Volume at 0.127mm Thick (mm^3)
Simple	681.06193	0.26813	0.13407
Tangential	485.74568	0.19124	0.09562
Meandering	631.43901	0.24860	0.12430
Fin Separator	474.58642	0.18685	0.09342
Flooded Flow	1225.17726	311.19502	155.59751

#### Flow Electrolysis Power Supply


Cat. No.	Description	Qty
F1030050	Flow Electrolysis Power Supply	Each

A full list of accessories and replacement parts can be found on our website



#### **Divided Cell (dual stream)**

In a divided cell the electrodes and their flow channels are separated by a PFSA Membrane, allowing for reactions to be performed that produce separate output streams. This can be advantageous if the two electrolyte solutions are incompatible or if the products of their reaction are troublesome to separate.

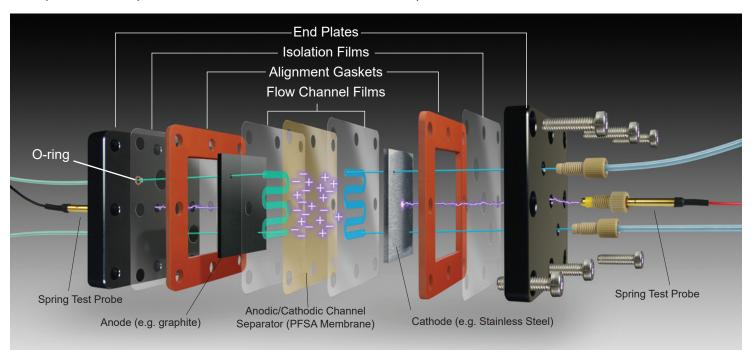




#### Flow Electrolysis Divided Cell (dual stream) Assembly

#### Assembly includes:

- Impervious Graphite Plate Electrode with holes (anode)
- 316L Stainless Steel Plate Electrode (cathode)
- · Flow Channel PFA Films (2 Meandering, 2 Tangential,
- 2 Simple, 2 Fin separator, 2 Flooded Flow)
- Isolation Layer PFA Films (2)
- PFSA Ion Exchange Channel Separator Membrane Orange Silicone Rubber Alignment Gaskets (2)
- Viton Chemical-resistant Internal O-rings, 2mm ID (4)
- Viton Chemical-resistant External O-rings, 3mm ID (4)
- Spring Test Probes/Pogo Pins (1 red, 1 black)
- Aluminum End Plates (2)
- Fittings and Hardware

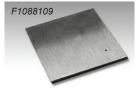



Cat. No.	Description	Qty
F1088200	Flow Electrolysis Divided Cell Ass. (dual stream)	Ea
Electro	de Plates	
F1088109	316L SS Plate Electrode (cathode)	Ea
F1088108	Impervious Graphite Plate Electrode (anode)	Ea
F1088208	Impervious Graphite Plate Electrode (anode) w/ Holes	Ea




#### **Consumable Accessories**

Cat. No.	Description	Qty
F1088405	PFA Isolation Layer Films	25
F1088504	Silicone Rubber Alignment Gaskets	10
F1088406	PFA Simple Channel Films	25
F1088403	PFA Meandering Channel Films	25
F1088407	PFA Tangential Mixer Channel Films	25
F1088410	PFA Fin Separator Channel Films	25
F1088412	PFA Flooded Flow Channel Films	25
F1088201	PFSA Ion Exchange Channel Separator Membrane	Each
F1088302	Viton Chemical-resistant External O-rings	50
F1088301	Viton Chemical-resistant Internal O-rings	50
F1088303	Spring Test Probes / Pogo Pins	10
F1030050	Flow Electrolysis Power Supply	Each





Accessories EquaVAP Evaporators Accessories

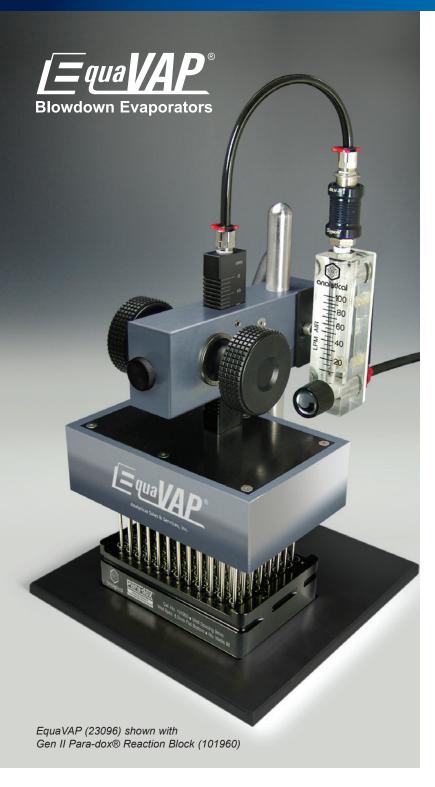

#### PRODUCT NOTE:

Requires regulated air supply set between  $80-110~\mathrm{psi}$ 

## EquaVAP® 96-Well Blowdown Evaporators - for Reaction Blocks

#### Rapidly evaporates common organic solvents

• Internal Flow Equalizers -


Distributes equal output across all needle ports to ensure even and symmetrical evaporation.

- Stainless Steel Needle Tips -Long term robust use, sturdy, wide solvent compatibilities
- Step-down Height Adjustment -Offers positive repeatable height positions
- Stand Alone -

Simple to operate, connects to Nitrogen or air supply with a standard socket

• Small Footprint (8" base) -

Fits into small glove boxes & standard fume hoods

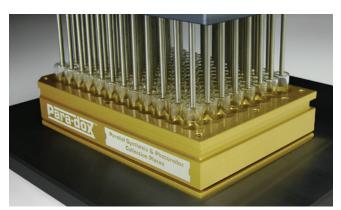

#### 96-Well EquaVAP Evaporator

 Use with 96-well Reaction Blocks with 1mL Vials or 96-well Polypropylene Collection Plates

С	at. No.	Description
2	3096	EquaVAP 96-Well Blowdown Evaporator


#### **EVAPORATION TIMES**

Chemical	Avg Evap Time* (HH:MM:SS)
Dichloromethane 99.9%	00:00:23
Acetone 100%	00:02:04
Methanol 99.9%	00:06:09
Ethyl Acetate	00:03:04
Acetonitrile	00:06:37
Isopropyl Alcohol 99.0%	00:08:48
Water (DI)	01:23:46

Conditions: 100 uL of solvent evaporated at AIR flow of 100 L/min (using 1mL round deep well plate)

\*Individual results may vary


#### EquaVAP® 48 and 24-Well Blowdown Evaporators - for Reaction Blocks

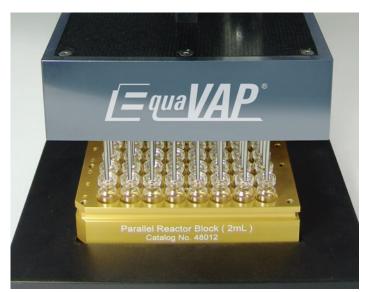

#### 48-Well EquaVAP Evaporator

 Use with 48-well Aluminum Reaction Blocks with 2mL Vials (101480), (48612), (48012)

Cat. No.	Description
23048	EquaVAP 48-Well Blowdown Evaporator for 48-Well Reaction
	Blocks with 2mL vials

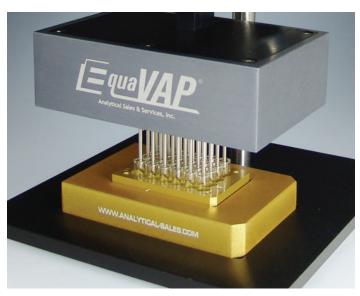

#### 24-Well EquaVAP for 9mm Vial Spacing

- Use with 24-well optimization blocks with 9mm vial spacing
- 24-well block fits into 96-well footprint with use of rack adapter (24245)

Cat. No.	Description
23024-09	EquaVAP 24-Well Evaporator for 24-Well Reaction Blocks with
	9mm Vial Spacing



#### 24-Well EquaVAP Evaporator (20mm)

- 20mm center-to-center needle distance
- Use with 24-well Aluminum Reaction Blocks with 1 Dram (4mL) or 2 Dram (8mL) Vials (24015), (24017), (24615), (24617)

Cat. No.	Description
23024-20	EquaVAP 24-Well Evaporator for 24-Well Reaction Blocks with
	20mm Spacing



#### 24-Well EquaVAP Evaporator (18mm)


- 18mm center-to-center needle distance
- Use with 24-well **Aluminum Reaction Blocks** with 18mm Well Spacing (101240), (24626)

Cat. No.	Description
23024-18	EquaVAP 24-Well Evaporator for 24-Well Reaction Blocks with
	18mm Spacing






Accessories Collection Plates & Cap Mats Accessories

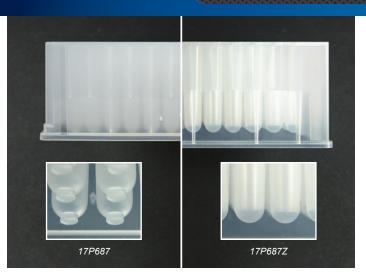

#### **Advantage™ 1mL Collection Plates**

#### Round Well with Round Bottom

Cat. No.	Description	Qty
17P687Z	1mL 96-Well Collection Plate with Round Well Bottoms	20
17P687	1mL 96-Well Collection Plate with Round Well Bottoms	20

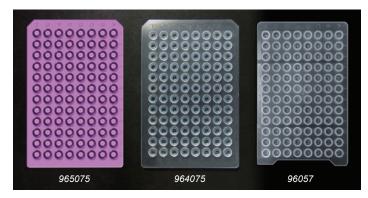

#### Advantage™ 96-Well Cap Mats - Round Well


Cat. No.	Description	Qty
965075	Purple Pre-Scored Ultra Thin Round Cap Mat	5
964075	Clear Pre-Slit Ultra Thin Round Well Cap Mat	5
96057	Autosampler Compatible 1mL Cap Mat with "X"	20



#### Advantage™ 2mL 96-Well Collection Plates Square Well with V-Shaped or Round Bottom

59623-23 2mL 96-Well Collection Plate with **V-Bottom** Wells

Cat. No.	Description	Qty
27P687	2mL 96-Well Collection Plate with Round-Bottom Wells	20

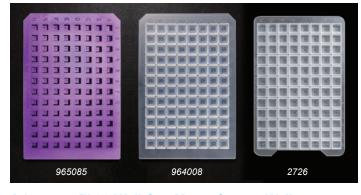

#### Advantage™ 96-Well Cap Mats - Square Well


Cat. No.	Description	Qty
965085	Purple Ultra Thin <i>Pre-Slit</i> Square Cap Mat	5
964008	Mighty Mat Purple Ultra Thin Square Cap Mat	5
2726	2mL Pierceable Cap Mat with "X" on Surface	20







## 24-Well Collection Plates and Cap Mats Cat. No. Description 24727 24-Well Collection Plate with Round-Bottom Wells. 10mL/well, 240m

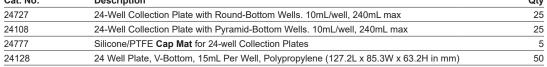



Fits nicely with Lumidox® II 24-well LED arrays








#### **Thermal Adapter Plates**

- For improved temperature transfer to samples
- For use with matching polypropylene collection plates

Our Thermal Adapter Plates conform to the well shape of their complementing polypropylene collection plate. They ensure uniform thermal transfer for heating and cooling when used in conjunction with a heating block or chiller (not included). The plate features a SLAS footprint and easy to identify purple anodized coating.

Analytical Sales & Services can make Thermal Adapter Plates to accommodate any shape and style polypropylene collection plate we offer. Please call 973-616-0700 for more information.

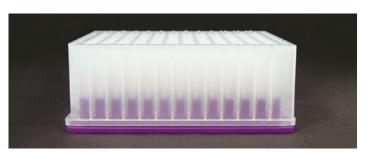



#### **Thermal Adapter Plates**

Cat. No.	Description	Qty
967720TAP	Thermal Adapter Plate for 967720 PP Collection Plate	Each
967720	2mL Round Well/Round Bottom 96-Well Collection Plate	25
59623-23TAP	Thermal Adapter Plate for 59623-23 PP Collection Plate	Each
59623-23	2mL Round Well/V-Bottom 96-Well Collection Plate	25
968810TAP	Thermal Adapter Plate for 1mL TrueTaper Collection Plate	Each
968810	1mL TrueTaper Collection Plate	25
968820TAP	Thermal Adapter Plate for 2mL TrueTaper Collection Plate	Each
968820	2mL TrueTaper Collection Plate	25
17P687ZTAP	Thermal Adapter Plate for 17P687Z PP Collection Plate	Each
17P687Z	1mL 17P687Z PP Collection Plate	25





Accessories Filter Plates Vial Trays and Loaders Accessories


#### **Vacuum Manifold Filtration System**

#### for SPE Sample Preparation

- Sturdy, clear acrylic construction
- Adjustable vacuum control
- Easy access to filtrate
- Compatible with all standard filter bottom microplates
- Collect into any storage plate
- · Compatible with robotic handling

The Advantage™ Vacuum Manifold Filtration System allows you to collect into any storage plate (including low volume collection plates) and is compatible with robotic handling systems. It uses an integral flat gasket between the collection plate and the filtration plate which stays totally secure within the system. A spacer is supplied for use with low volume collection plates. The Vacuum Manifold is constructed of tough acrylic for optimum performance and visual accessibility.

Cat. No.	Description	Qty
96844	Vacuum Manifold Filtration System for	Each
	SPF Sample Filtration	

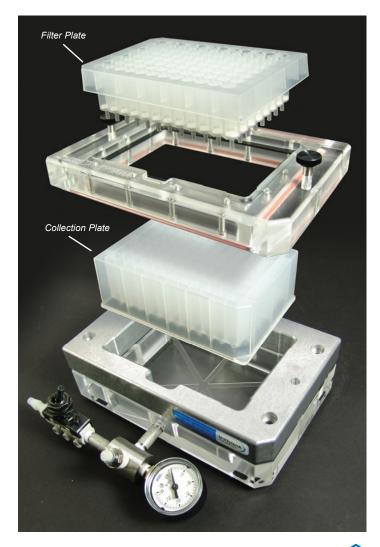



#### **Drug Discovery 96-Well Filter Plates**

• Low Protein Binding • General Filtration

Cat. No.	Description	Qty
964PP45	400μL Hydrophobic PP, 0.45μm	25

#### **General Filtration 96-well Filter Plates**

- After SPE Dilute and Shoot
- Use prior to LCMS and Microarraying

Low Protein Binding - Sample in Solvent

Cat. No.	Description	Qty
96245-10	2mL Hydrophilic PP, 0.45µm	10
96254-10	2mL Hydrophobic, PP 0.45μm	10

#### **Diatomaceous 96-Well Filter Plates**


- pH 1-13 No Pre-Treatment of the Bed is Necessary
- Packed with Flux Calcinated Diatomaceous Earth

Cat. No.	Description	Qty
96160-5	1mL Diatomaceous Filter Plate	5
96260-5	2mL Diatomaceous Filter Plate	5



**General Filtration 384-well Filter Plates** 

Cat. No.	Description	Qty
38407	384-Well Filter Plates, 140μL, Glass Fiber, 0.7μm	10
384603	Poly DVB SPE, 384 Well filter plate, 3mg/well	Each

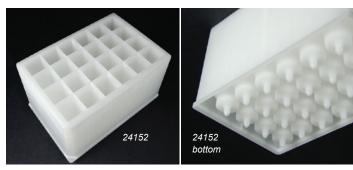

#### 24-well Filter Plates

Cat. No.	Description	Qty
24152	24 Well Filter Plates, PE 25µm, 15mL	15
2415SCX500	24 well SCX (Strong Cation exchange) 15 mL/500 mg bed	Each
241010	24 well Filter plates, Hydrophobic PP, 10 μm	25

#### **Assembled Vials in Stackable Trays**

- For easy loading Saves Time!
- Our thorough QC process ensures *Less Evaporation* compared to loose vials

#### 8x30mm Crimp Top Vials in Loaders







Recommended for top performance!

#### Shell Vials in Loaders



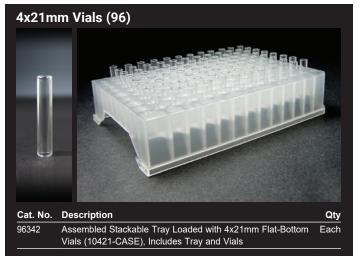

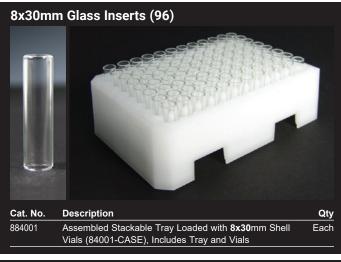





**Accessories** 

#### **Powder Transfer Plates**

Use to easily transfer powder from vial to vial or to mix powders between vials

- Wide funneled holes guide powder into vials
- · Grooves on bottom surface "lock" vials into position to avoid spillage





**Bottom Surface** 



## Powder Transfer Plate for 4x21mm Vials

#### Easily dispense powder into 4x21mm (50µL) Micro Vials

- 96-Position Polypropylene Tray with 4x21mm Vials (96342)
- 4x21mm Vials (10421-Case) in 96-Position Micro Optimization Block (96970)

**Powder** 

**Vials in Optimization Block (or Vial Holder Tray)** 

1. Lower plate onto vials and let the vials "settle" into the

grooves on the bottom surface of transfer plate

2. Powder can now be easily dispensed into each vial

Cat. No.	Description	Qty
96T100	Powder Transfer Plate for 4x21 Vials	Each

## Powder Transfer Plate with 4x21 Vials

#### Powder Transfer Plate for 8x30mm, 1mL Vials

Cat. No.	Description	Qty
96T105	Powder Transfer Plate for 8x30mm, 1mL Vials	Ea
884001	96-Position Stackable Tray with 8x30mm Shell Vials	Ea
	96T105	





with 8x30mm Vials (884001)

Use with 96-Position Stackable Tray loaded Flip repeatedly to mix powders between vials

#### Powder Transfer Plates for 15mm x 45mm, 4mL Vials

Cat. No.	Description	Qty
24T101	Powder Transfer Plate, 24-Well (with 20mm Spacing) for 15 x 45mm, 4mL Vials	Each
24T100	Powder Transfer Plate, 24-Well (with 18mm Spacing) for 15 x 45mm, 4mL Vials	Each



#### We are here to serve your needs quickly and efficiently! How do we do it?

- Personalized consultations and direct technical support
- Immediate response times
- Various convenient methods of contact
- Form submissions through our website
- Direct email (sales@analytical-sales.com)
- · Call us, a real person will answer!
- Virtual meetings with our sales team
  - Same day availability
  - Flexible scheduling
  - Accommodating international time zones
- Exceptional customer service
  - Receive a quote within less than one day
  - Email responses within hours, often minutes (during normal hours of operation)

#### **How to Order**

To place your order, you will need to supply the catalog number, a brief description and size, or the particular specifications when indicated.

#### You can order by:

Phone: 973-616-0700 • Fax: 973-616-0133

 Email: orders@analytical-sales.com Online store: www.analytical-sales.com

#### **Terms of Payment**

If you have an account with Analytical Sales, we will bill you for your purchases. All prices are F.O.B. Flanders, NJ. Terms of payment are net 30 days. To open an account, please call us. We also accept Visa®, Mastercard®, and American Express®. PayPal® is also available when ordering from our website.

#### Shipping

All items will be shipped via FedEx® Ground or common carrier unless otherwise instructed. Please examine all items immediately upon receipt. If you notice that an item was damaged in transit, it's important that you get a "damage notation" from the driver. If you notice damage upon unpacking an item, be sure to save all containers and packing materials. Please notify us immediately for instructions.

#### Returns

To receive credit for any product you return, you must first receive authorization. Please contact us for instructions. Returns must be made within 10 days of receiving authorization.

To see pricing please visit www.analytical-sales.com/catalog/pricing. All prices are subject to change without notice.







179 Rt 206 • Flanders, NJ • 07836

Phone: 973-616-0700 • Fax: 973-616-0133

Email: info@analytical-sales.com
Website: www.analytical-sales.com

